

PHD 2000 Syringe

Pump Series

User's Manual

<i>PHD 2000 Infusion</i>	<i>70-2000</i>
<i>PHD 2000 Infusion/Withdraw</i>	<i>70-2001</i>
<i>PHD 2000 Programmable</i>	<i>70-2002</i>
<i>PHD 2000 VHP</i>	<i>70-2023</i>



HARVARD

A P P A R A T U S

Table of Contents

SUBJECT	PAGE
1. Warranty and Repair Information	4
Manual Description	4
Warranty	4
Repair Facilities and Parts	4
Serial Numbers	4
Calibrations	4
General Safety Summary	5
2. Specifications and Ordering Information	6-7
3. Theory of Operation	8
4. Features	9
Pressure and Speed	9
Built-in Syringe Table and Custom Syringes.....	9
Infusion and Refill Rates	9
Target Volume	9
Auto Fill (Programmable Mode Only)	9
Modes of Operation.....	9
External Connections	10
Nonvolatile Memory	10
Stall Detection	10
Visual / Audible Alarm	10
Power-Up Options (Pump Mode Only)	10
Program Storage (Programmable Mode Only)	10
5. User Interface	11
Description of Keys	12
Entering Data	13
Special Features Access Table	13

Table of Contents

SUBJECT	PAGE
6. Initial Setup	14
7. Operation	15-20
Syringe Loading	15
Running the Pump	18
Diameter	18
Infuse Rate	18
Refill Rate	19
Target Volume	19
Auto Fill	19
Selecting the Run Mode	20
Pump Mode.....	20
Volume Mode	20
8. Program Mode	21-24
Program Description	21
Entering a Program.....	21
Sequence Operation	21
Program Printout.....	24
Program Run Time Error Messages	24
9. External Control and Interfaces	25-26
RS-232 Devices	25
Configuring Pump for RS-232 Devices.....	25
Pump Chain	25
Scale	26
Printer	26
TTL Devices	26
Foot Switch or Relay.....	26
Timer	26
Pumping Direction.....	26
Valve Control.....	26

Table of Contents

SUBJECT	PAGE
---------	------

10. Programming Tutorial	27
Multiple Infusion Example	27
Ramping Up Infusion Rate Example	28
Multiple Dispensing Example	29
Periodic Dispense Loop Example	30
Comb. of Infusion and Withdraw Profiles Ex. ...	31
Use of Events	32
11. Pump Chain Commands	34-45
Model 22 Protocol	35
Model 44 Protocol	37
A Table of Syringe Diameters	46

APPENDICES	PAGE
------------	------

B Stainless Steel Syringes	47
C Nominal Min./Max. Flow Rates	48
D PHD 2000 to PC Connection	49
E TTL Connector Specifications	49
F RS-232 Specifications	50
G Maintenance and Troubleshooting	51
H Accessories and Spare Parts	52
I Kits and Upgrades	53
J Custom Applications	54
K Symphony	54

Warranty and Repair Information



CAUTION: Verify operating voltage selection before plugging in the PHD2000 Pump. Refer to Section 6, Initial Setup.

Manual Description

This manual is designed to provide all operational and program information required to operate and maintain all three models of the PHD 2000 series pumps. The three models are: Infusion, Infuse/Withdraw and Programmable. The functions and features of these models are described in the Technical Specifications section.

Warranty

Harvard Apparatus warranties this instrument for a period of one year from date of purchase. At its option, Harvard Apparatus will repair or replace the unit if it is found to be defective as to workmanship or material. This warranty does not extend to damage resulting from misuse, neglect or abuse, normal wear and tear, or accident. This warranty extends only to the original customer purchaser.

IN NO EVENT SHALL HARVARD APPARATUS BE LIABLE FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES. Some states do not allow exclusion or limitation of incidental or consequential damages so the above limitation or exclusion may not apply to you. **THERE ARE NO IMPLIED WARRANTIES OF MERCHANTABILITY, OR FITNESS FOR A PARTICULAR USE, OR OF ANY OTHER NATURE.** Some states do not allow this limitation on an implied warranty, so the above limitation may not apply to you.

If a defect arises within the one-year warranty period, promptly contact *Harvard Apparatus, Inc. 84 October Hill Road, Building 7, Holliston, Massachusetts 01746* using our toll free number 1-800-272-2775. Goods will not be accepted for return unless an RMA (returned materials authorization) number has been issued by our customer service department. The customer is responsible for shipping charges. Please allow a reasonable period of time for completion of repairs, replacement and return. If the unit is replaced, the replacement unit is covered only for the remainder of the original warranty period dating from the purchase of the original device.

This warranty gives you specific rights, and you may also have other rights which vary from state to state.

Repair Facilities and Parts

Harvard Apparatus stocks replacement and repair parts. When ordering, please describe parts as completely as possible, preferably using our part numbers. If practical, enclose a sample or drawing. We offer a complete reconditioning service.

Serial Numbers

All inquires concerning our product should refer to the serial number of the unit. Serial numbers are located on the rear of the chassis.

Calibrations

All electrical apparatus is calibrated at rated voltage and frequency.



CAUTION: Not for clinical use on human patients.

Warranty and Repair Information

GENERAL SAFETY SUMMARY

Please read the following safety precautions to ensure proper use of your syringe pump. To avoid potential hazards and product damage, use this product only as instructed in this manual.

To Prevent Hazard or Injury:

Use Proper Line Cord

Use only the specified line cord for this product and make sure line cord is certified for country of use.

Ground the Product

This product is grounded through the grounding conductor of the power cord. To avoid electric shock, the grounding conductor must be connected to earth ground. Before making any connections to the input or output terminals of the product, ensure that the product is properly grounded.

Make Proper Connections

Make sure all connections are made properly and securely.

Observe all Terminal Ratings

Review the operating manual to learn the ratings on all connections.

Use Proper Fuse

Use only specified fuses with product.

Avoid Exposed Circuitry

Do not touch any electronic circuitry inside of the product.

Do Not Operate with Suspected Failures

If damage is suspected on or to the product do not operate the product. Contact qualified service personnel to perform inspection.

Place Product in Proper Environment

Review the operating manual for guidelines for proper operating environments.

Observe all Warning Labels on Product

Read all labels on product to ensure proper usage.



CAUTION
Refer to Manual



*Protective
Ground
Terminal*

Specifications

Specifications

	PHD 2000 SYRINGE PUMP SERIES	PHD 2000 VHP REMOTE SYRINGE PUMP SERIES
Accuracy	± 1%	± 1%
Reproducibility	± 0.1%	± 0.1%
Syringes (Min./Max.)	.5 µl / 140 ml	Holds 20 to 200 ml stainless steel syringes
Flow Rate:		
Minimum	.0001 µl/hr (with 0.5 µl Syringe)	1.5 µl/hr
Maximum	220.82 ml/min (with 140 ml Syringe)	112 ml/min
Calibration	Automatic	Automatic
Display	2 line, 40 char. fluorescent	2 line, 40 char. fluorescent
Non-Volatile Memory	Stores all settings	Stores all settings
RS-232 Interface	Multiplexed dual bidirectional ports	Multiplexed dual bidirectional ports
RS-232 Connectors	RJ11-4 cond. phone plug	RJ11-4 cond. phone plug
TTL Connectors	9 pin D-Sub. connector	9 pin D-Sub. connector
Linear Force (Max)		
Standard	23 kg (50 lbs)	196 kg (433 lbs)
High Force	30 kg (66 lbs)	N/A
Drive Motor	1.8° Stepper Motor	1.8° Stepper Motor geared
Motor Drive Control	Microprocessor from 1/2 to 1/32 microstepping	Microprocessor from 1/2 to 1/32 microstepping
Motor Step per one Rev. of Lead Screw	From 800 to 12,800	From 8,000 to 128,000
Step Rate:		
Minimum	27.3 sec/step	27.3 sec/step
Maximum	416.7 µsec/step	416.7 µsec/step
Pusher Travel Rate		
Minimum	0.18 µm/min	0.09 µm/min
Maximum	190.676 mm/min	71.210 mm/min
Power	115/230 VAC: 50-60 Hz External voltage selection 65 W, 0.5 A fuse	115/230 VAC: 50-60 Hz External voltage selection 65 W, 0.5 A fuse
Voltage Range	95-130 VAC, 220-260 VAC	95-130 VAC, 220-260 VAC
Dimensions, HxWxD	15.9 x 22.8 x 27.9cm (6.3 x 9 x 11in)	9.5 x 27.9 x 22.9 cm (3.75 x 11 x 9 in)
Syringe Holder, HxWxD	N/A	22.9 x 43.2 x 30.5 cm (9 x 17 x 12 in)
Remote Cable Length	N/A	152 cm (60 in) approx.
Weight	4.5 kg (10 lb)	

Ordering Information

Harvard PHD 2000 Syringe Pump Series				
Pump Versions	Standard (29 lbs Force)		High Power (58 lbs Force)	
	Regular	Remote	Regular	Remote
PHD 2000 Infusion Only	70-2000	70-2100	71-2000	71-2100
PHD 2000 Infuse/Withdraw	70-2001	70-2101	71-2001	71-2101
PHD 2000 Programmable Infuse/Withdraw	70-2002	70-2102	71-2002	71-2102
PHD 2000 Infusion Only with 6/10 Multirack	70-2003	70-2103	71-2003	71-2103
PHD 2000 Infusion Only with 4 x 140 Multirack	70-2004	70-2104	71-2004	71-2104
PHD 2000 Infusion Only with Microliter Rack	70-2005	70-2105	71-2005	71-2105
PHD 2000 Infuse/Withdraw with 6/10 Multirack	70-2006	70-2106	71-2006	71-2106
PHD 2000 Infuse/Withdraw w/ 4 x 140 Multirack	70-2007	70-2107	71-2007	71-2107
PHD 2000 Infuse/Withdraw with Microliter Rack	70-2008	70-2108	71-2008	71-2108
PHD 2000 Programmable with 6/10 Multirack	70-2009	70-2109	71-2009	71-2109
PHD 2000 Programmable with 4 x 140 Multirack	70-2010	70-2110	71-2010	71-2110
PHD 2000 Programmable with Microliter Rack	70-2011	70-2111	71-2011	71-2111

Harvard PHD 2000 Push/Pull Syringe Pump Series	
Pump Versions	Standard (29 lbs Force)
	Regular
PHD 2000 Infusion/Withdraw	70-2020
PHD 2000 Programmable Infuse/Withdraw	70-2019

Harvard PHD 2000 HVP (High Volume and Pressure)	
Pump Versions	Standard (over 400 lbs Force)
	Regular
PHD 2000 HVP (High Volume and Pressure)	70-2023

Theory of Operation

All three models in the PHD 2000 series employ a microcontroller which controls a small step angle stepping motor that drives a lead screw and Pusher Block. Micro-stepping techniques are employed to further reduce the step angle, eliminating flow pulsation. A keypad is used for entry of operating data to the pump. Data can also be entered via an RS-232 connector located on the rear panel. The microcontroller, using the internal Syringe Look Up Table, calculates the cross-sectional area of the syringe selected and calibrates the flow rate and volume accumulation. The numerous features of the PHD 2000 series result from the use of microprocessor technology.

The PHD 2000 Infuse/Withdraw model is equipped with all the functions of the PHD 2000 Infusion model along with additional features to provide the Withdraw (Refill) function. The PHD 2000 Infuse/Withdraw model's additional features include two additional function keys on the keypad, REFILL RATE key and the INFUSE/REFILL key and the addition of syringe Retaining Brackets and Adjustable Stops for refill. Use of the REFILL RATE key, the INFUSE/REFILL key, Retaining Brackets and Adjustable Stops are all described in detail in section 7 of this manual.

The PHD 2000 Programmable model is equipped with all the functions of the PHD 2000 Infuse/Withdraw model along with additional features to provide Programmable functions. The PHD 2000 Programmable models additional features include two additional function keys on the keypad, the PROGRAM key and the AUTO FILL key. Use of the PROGRAM key and the AUTO FILL key are described in detail in section 7 of this manual.

Pressure and Speed

PHD 2000 can deliver up to 220.82 ml/minute with a single 140 ml syringe. Maximum pressure is dependent on syringe size. For example a Becton-Dickinson 'Plasticpak' 50/60cc syringe run at 5ml/minute delivers an average maximum pressure of 1.8kg per cm² (25psi).

Built-In Syringe Table and Custom Syringes

If a non-standard syringe is to be used, enter the inside diameter of the syringe in millimeters. If a standard syringe is to be used, use the pump's built-in syringe table. Syringes are arranged according to manufacturer and material, and then according to size. The pump will look up and use the diameter for the syringe you select. (See Appendix A for a listing of the standard syringes).

Infusion and Refill Rates

Specify independent rates for infusing and refilling. This allows a slow infusion rate then a fast refill. If Refill Rate is not set, the software defaults to the set Infuse Rate value.

Target Volume

Specify the volume that is to be infused or refilled. The pump will run at the rate specified until this volume has been delivered when in the Volume mode.

Auto Fill (Programmable Model Only)

Auto Fill automatically activates an externally attached solenoid (Refer to Appendix H for part number) and refills the syringe when it is empty. This permits infusions to be virtually independent of syringe capacity.

Modes of Operation

Pump

Runs continuously, in the infuse or refill directions, until stopped. Refill not applicable to the Infusion model.

Volume

Runs until a specified volume has been pumped or refilled. Refill not applicable to the Infusion model.

Program

Pump operates according to a specified sequence of instructions. Programmable model only. (Note: All modes interact with Auto Fill feature, but it is only available on the Programmable model)

Features

External Connections

TTL

Allows pump operations to be synchronized with external devices or by a person at a distance from the pump. Direction of pump travel can be set via a TTL pin. Also, a TTL pin is used to control an external valve for refilling. Additional TTL pins are available for general use. (See Section 9 and Appendix E).

RS-232

Multiple pumps can be 'daisy chained' together and remotely controlled from a computer or any device communicating via RS-232. A scale can be connected, enabling the pump to infuse by weight instead of by volume. (Section 9). A printer can be connected to record final volumes or weights whenever the pump stops. In addition the program entered for the program mode can be listed on a connected printer. Both a scale and a printer may be connected simultaneously. (See Section 9).

Nonvolatile Memory

All operational data entered into the pump from the keypad or from a computer will be stored, including the program. On power up, the display will blink until the pump receives its first command and all settings from when it was powered down will be recalled.

Stall Detection

An optical detector is used to verify expected movement of the motor. If the motor is prevented from turning due to jamming or excessive back pressure, the pump will stop and the display will read: PUMP STALLED.

Visual/Audible Alarm

After a power interruption the alarm can be selected to be a silent flashing display or an audible signal in addition to flashing display.

Power-Up Options (Pump Mode only)

Enables the user to select between a standby mode (waiting for a command to begin running) after power is interrupted or a restart mode that begins running the pump at the settings that were in process when the power was interrupted.

Program Storage (Programmable model only)

Programmable model can store up to 4 sets of 9 program sequences for later selection.

User Interface

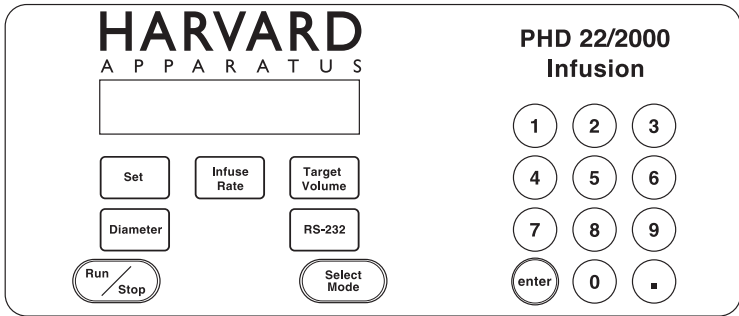


Figure 1. Infusion Model Keypad.

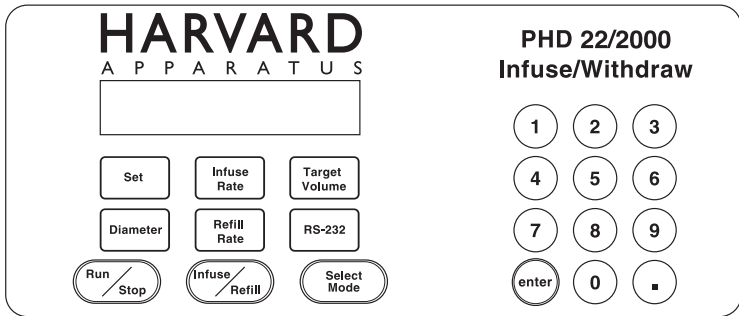


Figure 2. Infuse/Withdraw Model Keypad.

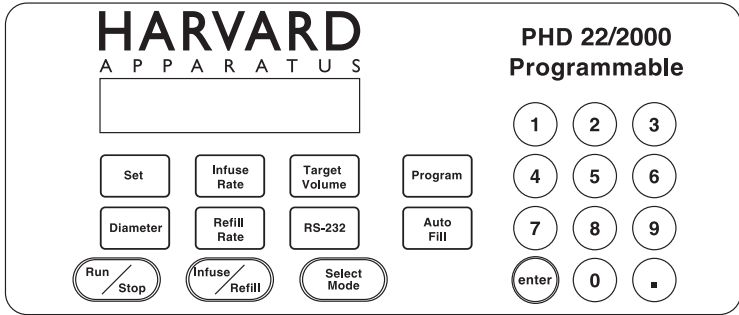


Figure 3. Programmable Model Keypad.

The user interface consists of a keypad with a display area. The display consists of a 2 line, 40 character alphanumeric vacuum fluorescent display. The display will show one of three types of messages: Default display, setting display, or an informational message. The top line of the display is divided into three sections. The first consists of the delivered volume, or time interval remaining in an operation, expressed with 5 digits plus a decimal point. Next is the units of the previous number expressed as 'ml' for milliliters, 'g' for grams, if a scale is attached, or 'time' for a time interval. Third is the pumping direction, either 'INFUSE' or 'REFILL', or the current operating state of the pump. A diamond symbol ◊ in the last character of the display indicates pump chain communication has been received.

User Interface

Setting displays are used to facilitate entering control information and data into the pump. Data entry will be discussed in detail later in this section. Informational messages occur at various times to indicate such items as a data setting out of range, or a detected problem, such as the pump stalling. Pressing any key clears the message from the display. The second line of the display is also divided into 3 sections. The first is the Running Mode (Pump, Vol, Prog). The second section is the AutoFill indication (no message if Auto Fill is Off, If On, AutoFill ON is displayed) The third section is a pair of directional arrow heads (< or >) indicating that the pump is running and in what direction. AutoFill is not applicable to the Infusion and Infuse/Withdraw Models.

DESCRIPTION OF KEYS

The keypad consists of 19 keys on the Infusion model, 21 keys on the Infuse/Withdraw model and 23 keys on the Programmable model. The keys are used for entering control information and data into the pump. These keys are grouped into 3 sections (see Figures 1, 2 & 3): Set keys, Toggle keys, and Data Entry keys.

Set Keys



SET – Allows modification of a data item in this group of keys. To modify a data item, press the relevant key after pressing the SET key. Pressing Set with an appropriate data entry key will also allow the selection of special features. Refer to Entering Data on page 7.

INFUSE RATE – Displays/sets current infuse rate. Scrolls through rate units with each press of the key when setting. Displays current programmed rate while running in program mode.

TARGET VOLUME – Displays/sets current volume mode target volume.

PROGRAM – Displays current program number, program sequence and sequence number. Held down, with a printer attached and the pump stopped, prints a program listing. Not applicable to the Infusion or Infuse/Withdraw models.

DIAMETER – Displays/sets current syringe diameter. When held down during setting, accesses built-in syringe table.

REFILL RATE – Displays/sets current refill rate. Scrolls through rate units with each press of the key when setting. Not applicable to the Infusion model.

RS-232 – Displays/sets current RS-232 device(s) attached.

AUTO FILL – Turns Auto Fill feature setting on/off. Also, displays/sets syringe refill volume.

Toggle Keys



Keys in this group, when permitted, successive states of the keys' function are selected when the key is pressed.

RUN/STOP – Starts/stops—interrupts running of pump.

INFUSE/REFILL – Changes direction of pusher block travel during operation in Pump and Volume modes. Pump must be stopped or in Pump Mode to reverse direction of pump. Not applicable to the Infusion model.

SELECT MODE – Toggles in an incremental loop through PUMP, VOLUME or PROGRAM run modes with each press of the key. The current mode is displayed on the 2nd line of the display. (Program mode is only available on the Programmable Model).

Data Entry Keys



1, 2, 3, 4, 5, 6, 7, 8, 9, 0, - – Used to enter numeric data values or access special features.

ENTER – Saves and stores displayed data value in memory when setting a data item.

User Interface

ENTERING DATA

Set Keys 

Keys with rectangular outlines in this group are used to modify or review settings of the pumps control data. To review the current setting of a control data item, simply hold down the relevant key and the data setting will appear in the top line of the display. In the case of the PROGRAM key (Programmable model only), if it is held down the data will be sent to the printer if one is attached.

To modify a data setting, first press then release the green outlined SET key. The display should then read 'SET WHAT?'. Press the key in the SET key group whose data is to be modified. The display will display the current setting. Data is entered into the pump by either entering a numerical value or by scrolling through a menu of choices. Always press the green outlined ENTER key to terminate and store each data request by the pump.

If you are to enter a numerical value, the far left of the display will show 'ENTER' followed by the units of the number to be entered. Using the numerical keys on the right side of the keypad (see Figures 1, 2 & 3) enter the new data value. Up to five digit numbers are accepted, including up to four decimal places. Entering more than five digits will clear the previous five digits entered in the display. Press the green outlined ENTER key when the desired data value is displayed. If the far left of the display does not show 'ENTER', then a menu of choices is being displayed. Pressing the relevant key, according to the choices being displayed, selects successive menu entries. When the desired selection is displayed, press the green outlined ENTER key.

If the data value entered is outside the pump's operating parameters, the display will read 'OUT OF RANGE'. Pressing any key will restore the display with the original data value. Enter another data value within the pump's parameters or just press the green outlined ENTER key to reuse the original data value.

The data value entered can be reviewed as described above. Note: Certain data items have multiple settings. For these, after the ENTER key is pressed, the display will prompt you for the additional information. Various rules apply to when, what and how data can be set at various times. See the relevant section for further details.

SPECIAL FEATURES ACCESS TABLE

Pressing the green outlined SET key and then the appropriate numerical key will allow the setting of several special features. Pressing the appropriate numerical key again will scroll between the options provided. Refer to the table below. When the desired option is displayed, press the green outlined ENTER key.

RS232 Protocol	SET	1
Alarm Mode	SET	2
Power Up Mode	SET	3

Initial Setup

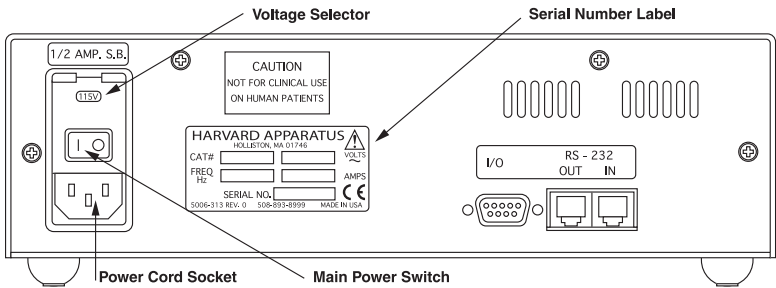


Figure 4. Rear Panel: Voltage Selector and Serial Number Label

1. Read the entire manual to become familiar with all features and functions of the PHD 2000.
2. If the operating voltage is other than 115VAC, 60Hz, locate the voltage selector on the rear panel (See Figure 4) of the pump and set it to the voltage being used. Also, if other than 115 V, 60 Hz is being used, the plug must be cut off and an appropriate plug installed, observing the polarity of the international line cord used:

Brown – live	Blue – neutral	Green – ground
--------------	----------------	----------------
3. Take this opportunity to fill out and mail the Warranty Card. The Catalog and Serial numbers are located on the Serial Number label on the rear panel. Refer to Figure 4.
4. Turn on main power switch (See Figure 4) located directly above the power cord receptacle on the rear panel. The two-line display will illuminate indicating that the power connections are correct. The flashing display indicates that power has just been applied. Pressing any key will reset the blinking display to constant illumination.
5. See Section 7 for operating Instructions on next page.
6. To change the Voltage setting from 115V to 230V, turn off the Main Power switch (0) and remove the power cord. Pry open the access door using a small screwdriver. Refer to (A) in Figure 4.
7. Remove the red fuse holder assembly by prying it out of the switch assembly using a small screw driver.
8. Rotate the red fuse holder 180° and install it back into the switch assembly. Close the small access door firmly until fully closed. 230V should be showing through the small opening in the switch assembly. Install the power cord.
9. Fuse Replacement. Turn off power, remove power cord from power module. Use straight blade screw driver, pry open access door. Remove fuse holder, remove fuse(s) from holder. Replace fuse(s) and fuse holder. (Replace holder with correct voltage marking is showing through access door window) For continued fire protection, replace fuse only with 250V fuse of the specified type and rating. (3AG 1/2 AMP SLO-BLO)



WARNING: Use in manner not specified by the manufacturer may impair the protection provided by the equipment.

Operation

Because of the wide range of functions that the PHD 2000 is capable of performing, certain information about your application must be entered into the pump. At minimum, the pump needs to know the Diameter of your syringe, the Infusion Rate and Direction of travel. (Direction of travel Infuse/Refill, is not applicable to the Infusion model) This is the only information needed to run the pump in the Pump Mode. If not specified, the Refill Rate (Refill Rate is not applicable to the Infusion model) will default to the Infuse Rate. The pump will need additional information to utilize its more advanced features. See Section 5, User Interface, for general information on data entry.


SYRINGE LOADING

Standard with a 2 Syringe Holder

The Pusher Block (2) and Syringe Holder Block (3) are fitted with movable retaining brackets (4 and 5) (Available as an option for the PHD 2000 Infusion model as an Anti-siphon kit - Catalog # 70-2015) which hold the syringe barrel and plunger firmly when refilling. When loading the syringe into the pump it is necessary to adjust these retaining brackets and the Adjustable Stops (8 & 9). The Pusher Block (2) is fitted with Release Tabs (1) to release the drive nut from the leadscrew so that the Pusher Block (2) can be moved freely on the leadscrew to facilitate loading the syringe. Refer to Figure 5.

1. Squeeze the Release Tabs (1) towards each other and move the Pusher Block (2) to the left allowing enough space for the syringe being used to fit between the Pusher Block (2) and the Syringe Holder Block (3).
2. Loosen the four screws to free the retaining brackets (4 and 5) on the Pusher Block (2) and Syringe Holder Block (3). This step is not required for the PHD 2000 Infusion model which is not equipped with retaining brackets (4 & 5) or Adjustable Stops (9).
3. Locate Syringe Clamp (6) and determine if clamp is in the correct position to firmly hold the desired syringe. Rule of thumb is that if the syringe is 20ml or larger, the tips of the clamp should be pointing down. If it is required to change the orientation of the clamp, unscrew the plastic knob located in the center of the clamp by turning it counter-clockwise (to the left). Remove the knob, lift the clamp off the shaft, flip the clamp over and place back on the shaft, then tighten the knob by turning it clockwise (to the right). Refer to Figure 6.
4. Lift up on the spring loaded Syringe Clamp (6) and rotate it 90° to allow clearance for the syringe. Place the barrel of the syringe on to the Syringe Holder Block (3) and rotate the clamp over the syringe. Ensure that the syringe barrel flange and plunger flange are located within the retaining brackets on both the Pusher Block (2) and the Syringe Holder Block (3). This step is not required for the PHD 2000 Infusion model which is not equipped with retaining brackets or Adjustable Stops (9). Lower the Syringe Clamp (6) down onto the syringe. CAUTION: Do not let the Syringe Clamp snap down on the syringe. Doing so may crack or break a glass syringe.
5. Ensure that both retaining brackets (4 and 5) are tight against the syringe barrel flange and plunger flange, then tighten the screws on the retaining brackets on the Pusher Block (2) and the Syringe Holder Block (3). This step is not required for the PHD 2000 Infusion model which is not equipped with retaining brackets or Adjustable Stops.
6. Squeeze the Release Tabs (1) and move the Pusher Block (2) to the right until the syringe plunger is located as far into the syringe as it will travel. Then squeeze the Release Tabs (1) and move the Pusher Block (2) to the left just enough so that the end of the syringe plunger is not touching the inside end of the syringe barrel and release the Release Tabs (1). This step is not required for the PHD 2000 Infusion model which is not equipped with retaining brackets or Adjustable Stops.

Operation

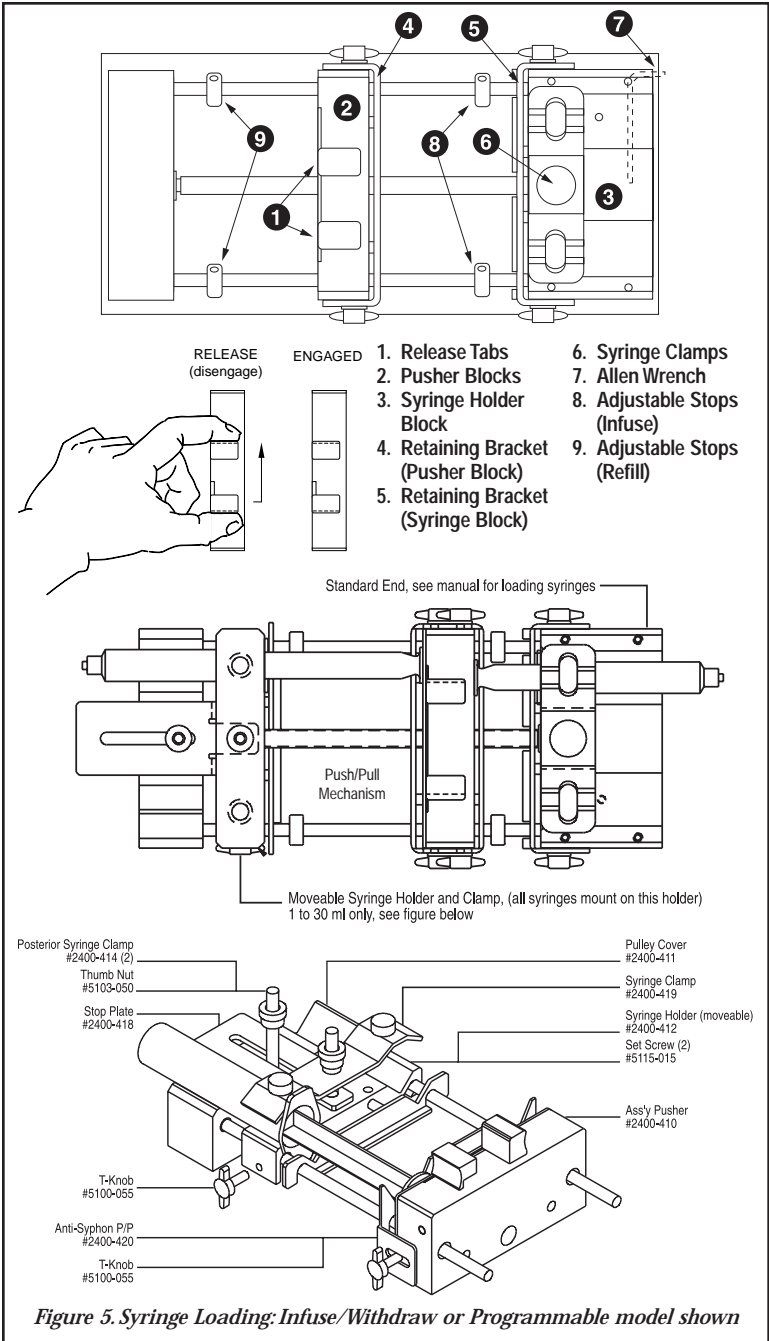

Operation

7. Locate Allen Wrench (7) in its holder on the back of the Syringe Holder Block and remove. Loosen front and rear Infusion Adjustable Stops (8) (right side) and slide them both to the left until they contact the retaining bracket (4). Using the Allen Wrench, tighten both Infusion Adjustable Stops (8).
8. Squeeze the Release Tabs (1) and move the Pusher Block (2) to the left until the syringe plunger is located as far out of the syringe as intended to travel. Then squeeze the Release Tabs (1) and move the Pusher Block (2) to the right just enough so that the end of the syringe plunger is just short of the maximum travel allowable and release the Release Tabs (1). This step is not required for the PHD 2000 Infusion model which is not equipped with retaining brackets or Adjustable Stops.
9. Loosen the front and rear Refill Adjustable Stops (9) (on the left side) and slide them both to the right until they contact the Pusher Block (2). Then using the Allen Wrench tighten both Refill Adjustable Stops (9) and return the Allen Wrench to its holder.

With Multiple Syringe Holder

Catalog Numbers - 70-2003, 70-2006, 70-2009

These Versions of the PHD 2000 have a saw tooth syringe holder capable of holding up to ten syringes, emptying them simultaneously at identical rates. This multiple syringe pump will accommodate up to:

- Ten syringes of any size up to 10 ml
- Six syringes of 20ml- 100 ml

Catalog Numbers - 70-2004, 70-2007, 70-2010

These Versions of the PHD 2000 have a multiple syringe holder which accommodates:

- Up to four syringes of 30ml to 140 ml

To load syringes rotate the spring loaded retainer arm out of the way and lay syringes in 'V' shaped slots with the syringe flange touching the holder. Rotate the retainer arm to hold syringes in place. If less than the maximum number of syringes are used, try to place them symmetrically about the lead screw to present a balanced load.

Maximum Force: The PHD 2000 will generate 29 lbs of force. With a 50 ml syringe it will deliver 25 psi.

Note: The back pressure for each syringe will be divided by the number of syringes used. Thus for four 50 ml syringes the maximum pressure developed in each syringe will be 6.3 psi.

Syringe Loading: Push/Pull Mechanism

1. Loading syringes on the standard end, see manual.
2. Loading syringes on push/pull end: 1 ml through 30 ml only.
Remove (2) thumb nuts and syringe clamp. Place syringes on moveable syringe holder. Adjust syringe retainer using (2) set screws. Position moveable holder for correct stroke. Clamp on to guide rod with the T-knob. Lock stop plate in place with thumb nut. Place syringe clamp over syringes and lock in place with thumb nut.

Operation

Running the Pump

Pressing the RUN/STOP key starts the pump. The pump will operate according to the relevant data entered as interpreted by the selected run mode. Pressing the RUN/STOP key while the pump is running stops the pump and the right side of the top line of the display will indicate 'INTERRUPT' plus '>' for infusing or '<' for refilling. (Refilling is not applicable to the Infusion model) This indicates that the pumping operation has been suspended and can be continued. Pressing the RUN/STOP key again will continue the pumping operation at the point that it was interrupted. Changing any of the settings, including the pumping direction and the run mode, cancels the interrupted operation and resets the display. The pump can also be started and stopped from a remote source.

Diameter

If the inside diameter of the syringe being used is known, enter the value in millimeters. Otherwise, access the built-in Syringe Table and select the syringe being used. After a new diameter is entered, directly or via the built-in Syringe Table, the Infuse Rate and Refill Rate are set to 0 and the Auto Fill feature is turned off (Refill Rate is not applicable to the Infusion model), (Auto Fill is applicable to the Programmable model only). This is done for reasons of safety. The maximum diameter that can be entered is 50 mm.

To access the built-in Syringe Table, after pressing the SET key then the DIAMETER key, hold down the DIAMETER key for about one second. Once the Syringe Table has been accessed, each press of the key will scroll through the manufacturers. Find the manufacturer and material, if applicable, of your syringe. Press the ENTER key to enter your selection. Now, using the DIAMETER key again, scroll by pressing the DIAMETER key until the size of your syringe is displayed, in cc or μL , as indicated on the display. Pressing the ENTER key will select the size of the syringe and look up and store the diameter. The diameter will be displayed until the ENTER key is released. Thereafter, pressing the DIAMETER key will display the selected diameter. In addition, the syringe size selected becomes the default Refill Volume when the Auto Fill feature is turned on. (Auto Fill is applicable to the Programmable model only) See Appendix A for a listing of the built-in Syringe Table and their respective diameters.

Infuse Rate

The Infuse Rate is the rate of pumping while infusing in the Pump or Volume modes. Also, the Infuse Rate is used as a starting rate for the program model if one is not specified in the program, regardless of pumping direction. (Programmable model only.) When entering the Infuse Rate, the INFUSE RATE key is used to scroll through the allowable units of rate. The allowable units are: ml/mn, ml/hr, $\mu\text{l}/\text{mn}$, $\mu\text{l}/\text{hr}$. While running in the Pump or Volume modes, the Infuse Rate can be changed. If the new rate is valid, it will take effect when the ENTER key is pressed. The minimum and maximum rates permitted vary depending on the diameter of the syringe. If an 'OUT OF RANGE' message is displayed when entering a rate, try using a different syringe for your application.

Operation

Refill Rate

The Refill Rate is the rate of pumping while refilling in the Pump or Volume Modes or during Auto Fill. If the Refill Rate hasn't been set (rate is 0), the Refill Rate will default to the Infuse Rate. (Refill Rate is not applicable to the Infusion model), (Auto Fill is applicable to the Programmable model only)

When entering the Refill Rate, the REFILL RATE key is used to scroll through the allowable units of rate. The allowable units are: ml/mn, ml/hr, µl/mn, µl/hr. (Refill Rate is not applicable to the Infusion model). While running in the Pump or Volume modes, the Refill Rate can be changed. If the new rate is valid, it will take effect when the ENTER key is pressed. (Refill Rate is not applicable to the Infusion model). The minimum and maximum rates permitted vary depending on the diameter of the syringe. If an 'OUT OF RANGE' message is displayed when entering a rate, try using a different syringe for your application.

Target Volume

The Target Volume is the volume that you desire to deliver at the set Infuse or Refill Rate. The pump must be set for this Volume mode for this feature to work. When entering the Target Volume, enter the volume desired in mls (milliliters) and press the ENTER key.

Auto Fill (only applicable to programmable model)

Use AUTO FILL key to toggle between Auto Fill 'ON' and 'OFF'. If Auto Fill is set to 'ON', the pump will next request the volume of the syringe in milliliters. The volume of the syringe is used as the refill volume of the syringe. This volume can be set to a volume smaller than the syringe if desired.

When set to 'ON', the syringe is assumed to be empty. Auto Fill continuously monitors the volume of the syringe according to the volume pumped. When the pump determines that the syringe is empty, the operation in progress is suspended and Auto Fill is activated. The pumping direction is then reversed and the pump runs at the refill rate. During the Auto Fill operation, the display will indicate the volume of liquid in the syringe.

When the volume in the syringe reaches the set refill volume, Auto Fill will stop, and the previous operation of the pump will resume. Auto Fill continues to monitor the volume of the syringe. TTL direction output is toggled 'ON' during refill. Refill Rate defaults to Infuse Rate if not set.

Note: Auto Fill will only activate while infusing. (i.e., if the pump direction is set to Refill, the pump will not stop when the syringe is full.) Also, if the syringe plunger is manually moved, the pump will lose track of the true syringe volume.

Operation

Selecting the Run Mode

After entering any necessary operating data into the pump, select the pumping mode that will be used when the pump is operated. Pressing the SELECT MODE key advances the second line of the display to indicate the next run mode. Advance the run mode to the desired mode, either Pump Mode, Volume Mode or Program Mode until it is displayed. (Program Mode is applicable to the Programmable model only)

Pump Mode

The pump will continuously pump, infusing or refilling, until stopped. While running, the Infuse and Refill Rates can be changed. The new rate, for the relevant pumping direction, takes effect when the ENTER key is pressed. Also, the pumping direction can be changed by pressing the INFUSE/REFILL key. (Refill, Refill Rate and Infuse/Refill are not applicable to the Infusion model),

Volume Mode

The pump will run, infusing or refilling, until a specified target volume is pumped or refilled. The TARGET VOLUME key is used to enter the Volume Mode target volume. Used in conjunction with Auto Fill, the target volume can be greater than the volume of the syringe. While running, the Target Volume, Infuse and Refill Rates can be changed. The new rate, for the relevant pumping direction, takes effect when the ENTER key is pressed. Target Volume can be used to enter a Target weight in grams if a scale is attached via RS-232. (Refill is not applicable to the Infusion model), (Auto Fill is applicable to the Programmable model only)

Program Mode

(only applicable to programmable model)

In the Program Mode the pump can make complex dispenses including changes in rate and target volume. These complex dispenses are easily programmed from the keypad and are detailed in the Program Mode and the Programming Tutorial sections.

Program Description

A program is made up of a set of sequences. Each sequence being a set of operating instructions for the pump to follow. When the pump is started in the PROGRAM run mode, the pump will start at sequence 1 and execute the operating instructions in that sequence. When the pump has completed the instructions for a sequence, it will go to the next, or specified, sequence and execute the instructions in that sequence. The pump continues this process until it either has reached a 'STOP' operation, the pump is manually or remotely stopped, or the last sequence has been completed. 4 programs with up to 9 sequences each may be entered.

A sequence consists of a sequence number, indicating the order of the sequence; a mode, indicating what operation the sequence will be performing; and the actual data for the operation, such as rates and volumes. The necessary data specified for each sequence will depend on the strategy used. One of two strategies may be chosen for a sequence's target. Strategy 1 (Target Volume) pumps until a target volume is reached, while Strategy 2 (Time Interval) pumps until a target time interval has lapsed. When Strategy 1 is used, enter a time interval of 0:00:00, then you will be prompted for the target volume. See the Programming Tutorial for example programs.

Entering a Program

It is advisable to plan out your program prior to entering the program into the pump. Press the SET key then the PROGRAM key to begin entering a program.

The following is a list of possible data that can be requested when entering a program and instructions on entering the data.

Sequence Operation

Use the PROGRAM key to select the sequence's operation. Operations that can be selected are: Profile, Increment (Incr), Decrement (Decr), Dispense, Event, Go To, Pause, Pump, TTL Out, Restart, Stop. When the required operation is displayed press the ENTER key. Additional information may be requested.

Rate

Enter the rate, using the INFUSE RATE key to change units. Note: If the rate entered is invalid, an error message will not be given at the immediate time of entry. An 'OUT OF RANGE' error message will be given during the running of the program.

Delta Rate

Enter the rate of Increment or Decrement desired. The units of the rate cannot be specified. Units will be the same as the units of the current pumping rate at the time the sequence is executed.

Target Volume

Enter the required delivered target volume of the sequence. For increment and decrement sequences, the target volume is an incremental target. An incremental target is added to the delivered volume at the start of the sequence.

Program Mode

Time Interval

Enter the time duration of the sequence in the form: 'hours : minutes : seconds'. If sequence Strategy 1 (Target Volume) is used, enter 0:00:00 for the time target. The maximum time interval is 9:99:99.

Number of Repetitions

Enter the number of times the sequence is to be repeated. The repetition number can be from 1 to 99,999.

Pumping Direction

Each sequence that specifies a pumping operation, also specifies a pumping direction. Use the INFUSE/REFILL key to change the pumping direction (< or >). (Direction of travel Infuse/Refill, is not applicable to the Infusion model)

Pin Level

Select either HI or LOW for the logic level of the programmable output pin 4. Use the PROGRAM key to change the setting.

Go to Sequence Number

Enter the destination sequence to continue operation of the program. Valid sequence numbers are 1 to 10.

Profile

Runs at specified flow rate until target volume is pumped or a time interval has elapsed.

Travel direction is as specified.

Data Specified:

Strategy 1:	Flow rate	Strategy 2:	Flow rate
(Volume)	Time Interval = 00:00:00	(Time)	Time interval
	Target volume		Pumping direction
	Pumping direction		

Incr

Increments current rate by specified value and pumps until the target volume is pumped or a time interval has elapsed. Units of rate will be that of the current rate of the pump or the infusion rate's units, if first sequence.

Sequence is repeated the specified number of times. Travel direction is as specified.

Data Specified:

Strategy 1:	Delta flow rate	Strategy 2:	Delta flow rate
(Volume)	Time Interval = 00:00:00	(Time)	Time interval
	Volume increment		Number of repetitions
	Number of repetitions		Pumping direction
	Pumping direction		

Program Mode

Decr

Same as INCR except rate is decremented.

Dispense

Repeatedly dispense specified volume. Runs at specified flow rate until a volume is pumped or a time interval has elapsed, then pump will stop. If no time interval was specified (Strategy 1), the display will show 'TRIGGER' and the next dispense will begin after an external or keyboard run command. Otherwise, the sequence will pause for specified time interval. Sequence is repeated the specified number of times. Travel direction is as specified.

Data Specified:

Strategy 1:	Flow rate	Strategy 2:	Flow rate
(Volume)	Time Interval 00:00:00	(Time)	Target volume
	Target volume		Time interval
	Number of repetitions		Number of repetitions
	Pumping direction		Pumping direction

Event

Program Events – A program event is an external event defined as a high to low transition on TTL pin-9. Within a program, a one time event trigger can be set which watches for and acts upon the external event. The triggered event causes an immediate continuation of the program at the specified sequence and the operation of the pump will be according to this sequence.

Data Specified:

Go To sequence number

Go To

Causes the program to immediately continue operation at the sequence specified.

Data Specified:

Go To sequence number

Pause

Pump stops for specified time then continues with next programmed sequence. Current program rate set to 0, with no change in units.

Data Specified:

Time interval

Pump

Runs the pump continuously at the specified flow rate without any pumping target. This mode can provide a background flow rate while waiting for an external event to trigger a new sequence specified by the EVENT operation.

Data Specified:

Flow rate
Pumping direction

Program Mode

TTL Out

Programmable TTL Pin: TTL output Pin 4 can be set to a HIGH or LOW level from within a program.

Data Specified:

TTL pin level

Restart

Immediately restart program from the first sequence.

Data Specified:

None

Stop

Stops pump and the program terminates.

Data Specified:

None

Program Printout

If a printer is attached and the pump is stopped, a program listing can be obtained by pressing the PROGRAM key for about one second. 'PRINTING PROGRAM' will be displayed while data is being sent.

Program Run Time Error Messages

If while running a program an operation is requested that cannot be performed, the pump will stop and an error message will be displayed. Error messages will be displayed with the following format:

Program N SEQ n: message

Where 'N' is the Program number and 'n' is the sequence number when an error was detected, and 'message' is the indicated error as follows:

INFINITE LOOP

A GO TO sequence cannot specify the current sequence.

INVALID GO TO

The target of the GO TO specified an invalid sequence number.

RATE UNDERFLOW

A decrement sequence decremented a rate to less than or equal to 0.

RATE OVERFLOW

An increment sequence caused an arithmetic overflow.

OUT OF RANGE

Specified or calculated rate is beyond the pumps capabilities with the specified syringe.

VOL TGT ERROR

A sequence with a volume target cannot follow a sequence with a time target, unless the volume delivered is zero or the pump is stopped at the start of the sequence entering it, select Program Mode using the SELECT MODE key and press the RUN/STOP key.

External Control and Interfaces

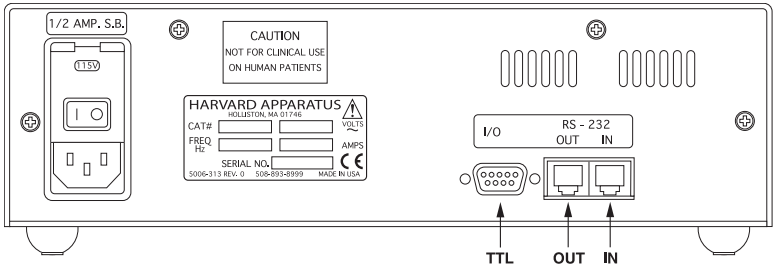


Figure 7. Rear Panel External Connections.

External devices that can be attached to the PHD 2000 are categorized into either RS-232 devices or TTL devices. Pump Chains, Scales and Printers are RS-232 devices, all other devices are TTL devices. See the Appendix F for specification details on attaching RS-232 devices and Appendix E for TTL.

RS-232 Devices

There are two telephone jack type connectors on the back of the unit (See Figure 7). These are the RS-232 ports. Looking at the back of the pump, the connector on the right is IN and the left is OUT. Attach the RS-232 connectors in the appropriate port according to the following chart:

Device	Port
Pump Chain computer side	IN
Pump Chain pump side	OUT
Scale	IN
Printer	OUT

Configuring the Pump for RS-232 Devices

Press the SET key and then press 1 to select either Model 22 or Model 44 Protocol (Refer to Section 11, Pump Chain Commands). Press the SET key, then use the RS-232 key to scroll through the menu of allowable RS-232 configurations. Possible configurations are:

- Pump Chain
- Scale
- Printer
- Scale & Printer

After entering the RS-232 configuration, additional information may be requested:

Pump Chain

Enter the 2-digit address assigned to the pump. Note: Each pump in the chain needs a unique address. After entering the address (up to 99), the baud rate will be requested. Use the RS-232 key to toggle between the supported baud rates: 1200, 2400, 9600 and 19200. Note: Each pump in the chain must have same baud rate. See the Section 11, Pump Chain Commands for pump chain control information.

External Control and Interfaces

Scale

Use the RS-232 key to toggle between the supported manufacturers: Mettler, Sartorius and Ohaus. When a scale is attached, the weight will be read from the scale and used as the delivered volume whenever the pumping direction of the pump is set to infuse. When refilling, the syringe diameter is used for volume calculations. When the scale weight is displayed, the units will be grams.

Printer

No additional information requested when entering. With a printer (capable of serial port communication) attached, the pump will print the delivered volume whenever the pump stops or the direction of pumping changes, except before and after Auto Fill of the syringe. If the pump stops due to the pump stalling, an asterisk (*) will be appended to the volume printed. In addition, the entered pump program can be listed on the printer by pressing the PROGRAM key for about one second, with the pump stopped.

TTL Devices

The pump does not need to be configured to attach a TTL device. Simply plug the device into the 9-pin connector on the rear of the pump. See Appendix E for wiring specifications.

Foot Switch or Relay

Used to start and stop the pump. Pressing the foot switch performs the same function as pressing the RUN/STOP key on the keyboard. The foot switch connector allows remote or automated operation of the pump.

Timer

Opening the timer input starts the pump. Closing the timer input stops the pump. The timer input allows for an externally controlled pumping interval.

Pumping Direction (Not applicable for Infusion model)

Sets the direction of pumping. Opening the directional input sets the pump to infuse. Closing the directional input sets the pump to refill. The pumping direction input is recognized only in the situations that the INFUSE/REFILL key would be recognized, i.e., when the pump is stopped or running in the Pump Mode.

Valve Control (Not applicable for Infusion Model)

The valve control output is an indicator of the direction of pump travel. When the output is high, the pump is set to Refill. A output low indicates Infuse.

Programming Tutorial

In the following examples, the diameter is 26.7 mm and the infuse rate is 50 ml/mn. To run a program after entering it, select Program mode using the SELECT MODE key and press the RUN/STOP key. The examples on the following pages include: Multiple Infusions, Ramping Up Infusion Rate, Multiple Dispensing, Periodic Dispense Loop, Combination Infuse and Withdraw Profile, Use of Events and Use of Signal.

Multiple Infusion Example

The following program will instruct the pump to infuse according to the above graph. The program instructs the pump to infuse 10 ml at 75.000 ml/mn then infuse another 5 ml at 25 ml/mn then stop, for a total of three sequences. Since this graph is Rate vs. Volume, Strategy 1 will be used when entering the program.

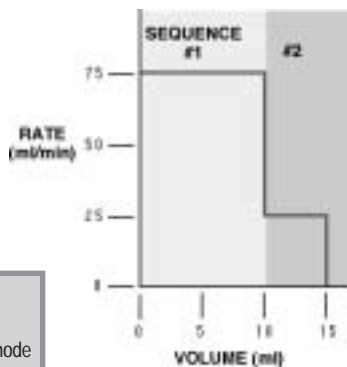


Figure 6. Multiple Infusions.

SEQUENCE 1:

Key Presses	Explanation
SET	Allows selection of Programming mode
PROGRAM	Press until PROFILE selected
ENTER	Enters selection
75	Enter rate of 75.000 ml/mn
INFUSE RATE	Press until units are ml/mn
ENTER	Enters rate
0 ENTER	Enter 0 for time, this indicates Strategy 1
10 ENTER	10 ml is the first target volume
INFUSE/REFILL	Toggles direction to infuse
ENTER	Enters sequence's pumping direction

SEQUENCE 2:

Key Presses	Explanation
PROGRAM	Press until PROFILE selected
ENTER	Enters selection
25	Enter rate of 25.000 ml/mn
INFUSE RATE	Press until units are ml/mn
ENTER	Enters rate
0 ENTER	Enter 0 for time, this indicates Strategy 1
5 ENTER	5 ml is the second target volume
INFUSE/REFILL	Toggles direction to infuse
ENTER	Enter sequence's pumping direction

SEQUENCE 3:

Key Presses	Explanation
PROGRAM	Press until STOP selected
ENTER	Enters selection and ends program entry

PROGRAM PRINTOUT

```
PROG1 SEQ 1: PROFILE
75.000 ml/mn
10.000 ml
INFUSE
PROG1 SEQ 2: PROFILE
25.000 ml/mn
5.0000 ml
INFUSE
PROG1 SEQ 3: STOP
```

Programming Tutorial

Ramping Up Infusion Rate Example

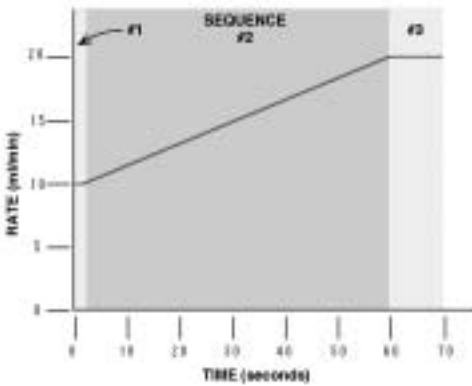


Figure 7. Ramping Up Infusion Rate

PROGRAM PRINTOUT

```

PROG1 SEQ 1: PROFILE
10.000 ml/mn
0:00:01 INTERVAL
INFUSE
PROG1 SEQ 2: INCR
0.1695 INCR
0:00:01 INTERVAL
INFUSE
59 REPEAT
PROG1 SEQ 3: PROFILE
20.000 ml/mn
0:00:10 INTERVAL
INFUSE
PROG1 SEQ 4: STOP
  
```

In this example, the pump will ramp up from 10 ml/mm to 20 ml/mn over 60 seconds, then continue to run at 20 ml/min for another 10 seconds. This is a Strategy 2 Program requiring four sequences:

1. Specify the initial rate as a profile of 10 ml/mn for one second.
2. Specify the ramp up to 20 ml/mn. Since the minimum resolution of an increment is one second, it will take 59 steps to reach the target rate. Sequence 2 starts at time 1 second and ends at time 60 seconds, giving it a duration of 59 seconds.

At one second a step, 59 seconds divided by one second per step equals 59 steps. The increase per step will be 20 ml/mn minus 10 ml/mn, divided by 59 steps or 0.1695 rounded to four decimal places.

3. Continue running at 20 ml/mn for 10 seconds with a profile operation.
4. Stop the pump.

Multiple Dispensing Example

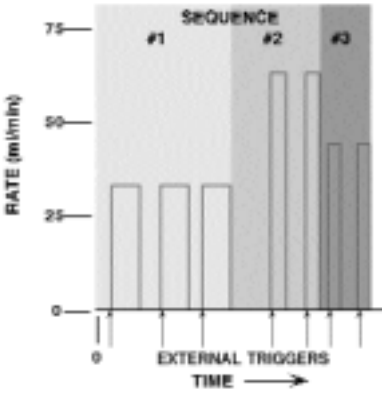


Figure 8. Multiple Dispensing

PROGRAM PRINTOUT

```
PROG1 SEQ 1: DISPENSE
35.000 ml/mn
15.000 ml
3. REPEAT
INFUSE
PROG1 SEQ 2: DISPENSE
65.000 ml/mn
25.000 ml
2. REPEAT
INFUSE
PROG1 SEQ 3: DISPENSE
45.000 ml/mn
17.000 ml
2. REPEAT
INFUSE
PROG1 SEQ 4: STOP
```

Here, a series of dispenses are programmed. Each dispense is started by a trigger, such as pressing the RUN/STOP key or pressing an attached foot switch. Seven dispenses are programmed: three of 15 ml at 35 ml/mn, two of 25 ml at 65 ml/mn, and two of 17 ml at 45 ml/mn. The pump's display will show 'TRIGGER' when it is waiting for a run trigger and < or > will not be displayed.

This is a Strategy 1 dispense. A time interval of 0 is specified when entering a Strategy 1 dispense. Since the total volume to be dispensed is 129 ml and the syringe volume is 50 ml, the Auto Fill feature would be very useful with this program.

Programming Tutorial

Periodic Dispense Loop Example

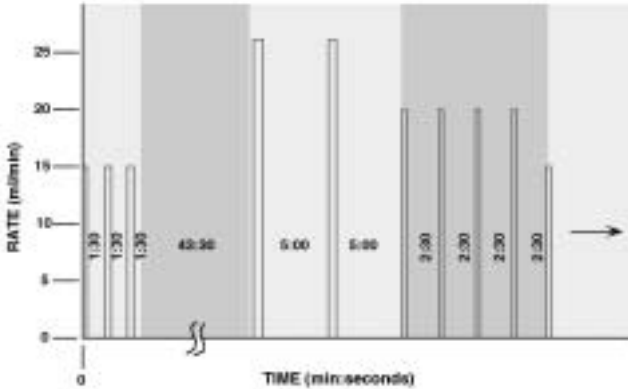


Figure 9. Periodic Dispense Loop

This is an example of a series of periodic dispenses of varying volumes and intervals. For this application, Strategy 2 dispenses are used. Note that between the third and fourth dispenses is a 45 minute interval. Each dispense in the first sequence is separated by a pause interval of 1:30. Since after the third dispense there already will be a 1:30 pause, an additional pause of 43:30 is used to extend the pause to the desired 45:00. Sequence 5 is a RESTART command, causing the series of dispenses to be continuously repeated until the pump is stopped.

PROGRAM PRINTOUT

```

PROGRAM PRINTOUT
PROG1 SEQ 1: DISPENSE
15.000 ml/mn
3.5000 ml
0:01:30 INTERVAL
3. REPEAT
INFUSE
PROG1 SEQ 2: PAUSE
0:43:30 INTERVAL
PROG1 SEQ 3: DISPENSE
25.700 ml/mn
6.7500 ml
0:05:00 INTERVAL
2. REPEAT
INFUSE
PROG1 SEQ 4: DIS-
PENSE
20.000 ml/mn
4.3000 ml
0:02:30 INTERVAL
4. REPEAT
INFUSE
PROG1 SEQ 5: RESTART
  
```

Combination of Infusion and Withdraw Profiles

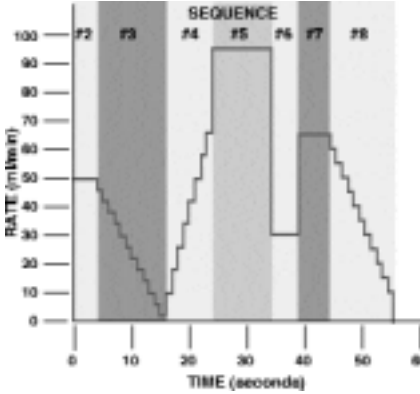


Figure 10. Combination Infusion and Withdraw Profiles

Example

Here is an example of a more complex profile program. Each 'run' of the infusion has been determined to pump 43.155 ml. The first sequence refills the syringe with the volume to be infused then the infusion profile is started, after which the syringe is refilled and the infusion is repeated until the pump is stopped.

PROGRAM PRINTOUT

PROG1 SEQ 1: PROFILE 75.000 ml/mn 43.155 ml REFILL PROG1 SEQ 2: PROFILE 50.000 ml/mn 0:00:04 INTERVAL INFUSE PROG1 SEQ 3: DECR 4.0000 DECR 0:00:01 INTERVAL 12. REPEAT INFUSE PROG1 SEQ 4: INCR 5.0000 DECR 8.0000 INCR 0:00:01 INTERVAL 8. REPEAT INFUSE	PROG1 SEQ 5: PROFILE 95.000 ml/mn 0:00:10 INTERVAL INFUSE PROG1 SEQ 6: PROFILE 30.000 ml/mn 0:00:05 INTERVAL INFUSE PROG1 SEQ 7: PROFILE 65.000 ml/mn 0:00:05 INTERVAL INFUSE PROG1 SEQ 8: DECR 5.0000 DECR 0:00:01 INTERVAL 11. REPEAT INFUSE PROG1 SEQ 9: RESTART
---	--

Use of Events

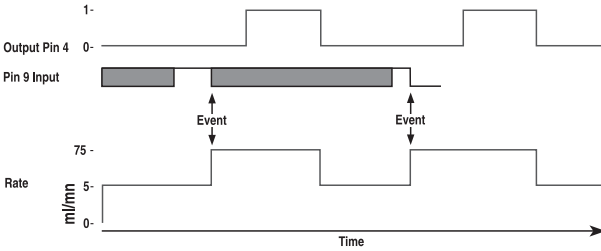


Figure 11. Use of Events

This is an example of the PHD 2000 working interactively with other laboratory equipment. The pump will continuously pump at 300 ml/hr until an external event, a high to low transition at pin 9, possibly produced by another PHD 2000, causes the pump to deliver a 15 ml bolus at 75 ml/mn. After delivering 5 ml of the bolus, output pin 4 is set to a logic high for the duration of the bolus after which it is dropped. This output pin can be attached to the timer input of another pump, such as a Harvard Pump 22 or 44, to create a precise mixture during the bolus. After the bolus is completed, the Pump 22 would be stopped and the PHD 2000 would return to delivering its background rate of 300 ml/hr, waiting for another external event.

PROGRAM PRINTOUT

PROG1 SEQ 1: TTL OUT OFF	5.0000 ml INFUSE
PROG1 SEQ 2: EVENT GO TO 4	PROG1 SEQ 5: TTL OUT ON
PROG1 SEQ 3: PUMP 300.00 ml/hr INFUSE	PROG1 SEQ 6: PROFILE 75.0000 ml/mn 10.0000 ml
PROG1 SEQ 4: PROFILE 75.0000 ml/mn	INFUSE PROG1 SEQ 7: RESTART

Programming Tutorial



Figure 12. Use of TTL Signal

Before the pumping flow begins, the pump sends a TTL signal to another piece of equipment for 5 seconds, but first verifies that the TTL line is low by turning it off for 1 second, then turns it on. The pumping process consists of an alternation of pumping 3 ml's at 53 ml/mn and 5 ml's at 75 ml/mn. It uses the TTL output to signal to another device whose sequence it is executing by lowering the TTL line before sequence 5 and raising it before sequence 8.

PROGRAM PRINTOUT

PROG1 SEQ 1: TTL OUT	53.000 ml/mn
OFF	3.0000 ml
PROG1 SEQ 2: PAUSE	INFUSE
0:00:01 INTERVAL	PROG1 SEQ 7: TTL OUT
PROG1SEQ 3: TTL OUT	ON
ON	PROG1 SEQ 8: PROFILE
PROG1 SEQ 4: PAUSE	75.000 ml/mn
0:00:01 INTERVAL	5.0000 ml
PROG1 SEQ 5: TTL OUT	INFUSE
OFF	PROG1SEQ 9: GO TO 5
PROG1SEQ 6: PROFILE	

This is an example of the PHD 2000 being operated from a remote location. When the pump is powered on, the position of the pusher block is unknown and must be homed to a known position.

At the syringe full position, a limit switch is placed such that it is tripped by the pusher block when the syringe is full. The limit switch is connected to pins 4 and 9, programmable output and the event input, on the TTL connector. When the pump receives a start signal, it first refills the syringe and stops when the limit switch is sensed. The pump then waits for a start trigger and performs a dispense then refills the syringe and waits again for the next start trigger.

PROGRAM PRINTOUT

PROG1 SEQ 1: EVENT	75.000 ml/mn
GO TO 7	PROG1 SEQ 7: DISPENSE
PROG1 SEQ 2: TTL OUT	10.000 ml/mn
ON	0.0001 ml
PROG1 SEQ 3: PAUSE	1. REPEAT
0:00:01 INTERVAL	REFILL
PROG1 SEQ 4: TTL OUT	PROG1 SEQ 8: PUMP
OFF	75.000 ml/mn
PROG1 SEQ 5: PAUSE	10.000 ml
00:00:01 INTERVAL	INFUSE
PROG1 SEQ 6: PUMP	PROG1 SEQ 9: RESTART

Pump Chain Commands

The Pump Chain RS-232 interface is used to enhance the control applications of the PHD 2000. This interface allows all control information, including a program, to be entered into the pump from an external source such as a computer. In addition, this interface allows up to 100 pumps and, in certain cases, other RS-232 devices to be controlled from a single RS-232 communication port on a computer.

Assign each pump in the pump chain a unique address from 0 to 99. This address is used to identify which pump is to receive a command and which pump is responding. Configure each pump with its assigned address and the baud rate being used (See Section 9, External Control & Interfaces).

When the control device communicates with the pump, a diamond appears on the far right of the display indicating that it is receiving RS-232 commands. The diamond remains on the display until the pump is turned off or SET RS-232 is entered on the keyboard, indicating a change in the RS-232 configuration.

A pump will not respond to pump chain communication while it is in a setting mode (entered when user presses the SET key). The pump can still be controlled from the keyboard while it is in a pump chain. Control data that is changed via RS-232 will be stored in the pump's non-volatile memory.

After each command is received and executed, the pump terminates its responses with a prompt. A prompt is a string of ASCII characters sent by a pump.

The PHD 2000 is capable of using two sets of pump interface commands. This will allow users of Harvard Apparatus Model 22 and Model 44 pumps to use existing programs to control the PHD 2000 pumps. The two sets of commands are contained on the following pages.

Pump Chain Commands

Each command sent to the pump chain is a string of ASCII characters, with leading zero's on numbers and all spaces optional. Numbers are a maximum of five digits. The following symbols are used in describing the commands:

Pump Chain Commands - Model '22' Protocol (Model 44 Protocol on page 37)

After each transmission to the Pump terminating with a CR character (ASCII 13), the pump enters remote mode and responds with the three character sequence:

CR LF prompt

The prompt character indicates the status of the pump as follows:

:	When stopped	(ASCII 58 decimal)
>	When running forward	(ASCII 62 decimal)
<	When running reverse	(ASCII 60 decimal)
*	When stalled	(ASCII 42 decimal)

Note: With a pump chain, if no address precedes the command transmitted, the pump at address 0 will respond to the command.

Commands - Response: CR LF prompt

RUN	Infuse (forward direction)
REV	Start (reverse direction) Not accessible on Infusion model
STP	Stop
CLV	Clear volume accumulator to zero
CLT	Clear target volume to zero
MLM number	Set rate, units are milliliters per minute
ULM number	Set rate, units are microliters per minute
MLH number	Set rate, units are milliliters per hour
ULH number	Set rate, units are microliters per hour
MMD number	Set diameter, units are mm. Rate is set to 0
MLT number	Set target infusion volume, units are ml.

Numbers between 0 and 1999 will be accepted by the pump. Leading zeros and trailing decimal point are optional. Any number of digits to the right of the decimal point may be transmitted. The number received will be rounded to four significant digits if the leading digit is 1, or three significant digits if the leading digit is 2 to 9.

Pump Chain Commands

Queries - Response: CR LF Value CR LF Prompt

Queries with Numeric Response

DIA	Send diameter value, units in mm
RAT	Send rate value in current range units
VOL	Send current accumulated infused volume, units are ml.
TAR	Send target volume, units are ml.
VER	Send model and version number

Value format: nnnn.nnn

The transmitted value is an 8 character string with leading zeros converted to SP characters. (ASCII 32) The fifth character is a decimal point (ASCII 46).

Query - Response: CR LF Range CR LF Prompt

Queries with String Response

RNG	Send range message
-----	--------------------

Range is a character string, one of: ML/H ML/M UL/H or UL/M

Error Responses

CR LF ? CR LF prompt	Unrecognized command
CR LF O O R CR LF prompt	Out of range

Pump Chain Commands

Pump Chain Commands - Model '44' Protocol

<u>Symbol</u>	<u>Meaning</u>
[...]	optional
{...}	select one
	either/or
f	digits 0 – 9 or a decimal point
d	digits 0 – 9
<cr>	carriage return (ASCII 13)
<lf>	line feed (ASCII 10)
<float>	f f f f f
<integer>	dddd
<time>	d:dd:dd
<text>	any string of ASCII characters

Command Formats and Meanings:

<u>Command Format</u>	<u>Meaning</u>
<cr>	Stops all pumps. All pumps on the pump chain interpret this as a stop command.
pump address, <cr>	Request for prompt The pump with the indicated address responds with its prompt
optional pump address, command, <cr>	Send a command to a pump. The pump with the indicated address executes the command then responds with its prompt. The optional pump address, if not specified, will default to pump address 0.

After each command is received and executed, the pump acknowledges the command with a prompt. Preceding the prompt may be some additional text responses. The additional text will be one or more lines of ASCII text, each preceded by a line feed and terminated by a carriage return:

<lf>, <text>, <cr>

A prompt is a string of ASCII characters sent by a pump indicating the pumps address and its present state:

<lf>, 1 or 2 digit address, prompt character

<u>Prompt Characters</u>	<u>Meaning</u>
:	Pump stopped
>	Pump infusing
<	Pump refilling
/	Pause interval (pump stopped)
*	Pumping interrupted (pump stopped)
^	Dispense trigger wait (pump stopped)

Pump Chain Commands

Pump Commands and Responses

- RUN** Starts pumping according to the present setting of the pump. If pump is already pumping, a 'Not Applicable' response will be given.
- STP** Stops pump if it was running. If pump was already stopped, a 'Not Applicable' response will be given.
- DEL** Request for volume delivered, in ml.
Response is of the following format:
space, space, f f f f f
- CLD** Request to zero volume delivered. If the pump was interrupted, it will cancel the interrupted condition. If the pump is running, request will not be accepted and a 'Not Applicable' response will be given. Otherwise, no response is given.

RAT [<float> [<units>]]

Request to set or query infusion rate setting.

Set infusion rate:
RAT rate

Set infusion rate and units:
RAT rate units

Rate is of format: f f f f f

Units are one of:

	<u>Definition</u>
UM	µl/mn
UH	µl/hr
MM	ml/mn
MH	ml/hr

If rate is accepted and valid, rate will become the new infusion rate.

If the rate is invalid, an 'Out Of Range' response will be given.

Command will not be accepted if the pump is running in the Program Mode and

a 'Not Applicable' response will be given.

Query infusion rate: RAT

Response is of the following format:
space, space, f f f f f units

Where units is one of the following:

ml/mn
ul/mn
ml/hr
ul/hr

Pump Chain Commands

RFR [<float> [<units>]]

Request to set or query refill rate setting.

Set refill rate:

RFR rate

Set refill rate and units:

RFR rate units

Rate is of format: f f f f f f

Units are one of:

Definition

UM	µl/mn
UH	µl/hr
MM	ml/mn
MH	ml/hr

If rate is accepted and valid, rate will become the new refill rate.

If the rate is invalid, an 'Out Of Range' response will be given.

Command will not be accepted if the pump is running in the

Program Mode and

a 'Not Applicable' response will be given.

Query refill rate: RFR

Response is of the following format:

space, space, f f f f f f units

Where units is one of the following:

ml/mn
ul/mn
ml/hr
ul/hr

PGR Request for the rate of pumping set during the running of a program.

Response is of the following format:

space, space, f f f f f f units

Where units is one of the following:

ml/mn
ul/mn
ml/hr
ul/hr

Pump Chain Commands

DIA [<float>]

Request to set or query syringe diameter setting.

Set diameter:

DIA diameter

Diameter is of format: f f f f f f

Units are MM.

INFUSE and REFILL rates will be set to zero and AUTO FILL will be set to off.

If diameter is accepted and valid, diameter becomes new diameter.

Diameter will not be accepted if the pump is running and a 'Not Applicable' response will be given.

If the diameter is invalid, an 'Out Of Range' response will be given.

Query diameter: DIA

Response is of the following format:

space, space, f f f f f f

Units are MM.

TGT [<float>]

Request to set or query target volume setting.

Set target volume:

TGT volume

Volume is of format: f f f f f f

Units are ML.

If volume is accepted and valid, volume becomes new target volume.

Volume will not be accepted if the pump is running and a 'Not Applicable' response will be given.

If the volume is invalid, an 'Out Of Range' response will be given.

Query volume: TGT

Response is of the following format:

space, space, f f f f f f

Units are ML.

Pump Chain Commands

MOD [(PMP | VOL | PGM)]

Request to set or query pumping mode

Set: MOD PMP (Puts pump in Pump Mode)
 MOD VOL (Puts pump in Volume Mode)
 MOD PGM (Puts pump in Program Mode)

Command will not be accepted if the pump is running and a 'Not Applicable' response will be given.

Query: MOD

If mode is PUMP, response will be:

PUMP

If mode is VOLUME, response will be:

VOLUME

If mode is PROGRAM response will be:

PROGRAM

DIR [(INF | REF | REV)]

Request to set or query pumping direction

Set: DIR INF (sets pumping direction to infusion)
 DIR REF (sets pumping direction to refill)
 DIR REV (reverses current pumping direction)

Command will not be accepted if the pump is running in volume or program modes and a 'Not Applicable' response will be given.

Query: DIR

If pump direction is infusion, response will be:

INFUSE

If pump direction is refill, response will be:

REFILL

AF [(ON | OFF)]

Request to set or query auto fill setting

Set: AF ON (turns Auto Fill feature on)
 Note: The syringe volume is also needed for auto fill to operate.
 (See SYR command)

AF OFF (turns Auto Fill function off)

Command will not be accepted if the pump is running and a 'Not Applicable' response will be given.

Query: AF

If Auto Fill function is ON, response will be: ON

If Auto Fill function if OFF, response will be: OFF

Pump Chain Commands

42

SYR [<float>]

Request to set or query syringe volume setting for auto fill.
Used in conjunction with Auto Fill feature. (See AF command).

Set syringe volume: SYR volume

Volume is of format: f f f f f

Units are ML.

If volume is accepted and valid, volume will become the new syringe Auto Fill volume.

Volume will not be accepted if the pump is running and a 'Not Applicable' response will be given.

If the volume is invalid, an 'Out Of Range' response will be given.

Query syringe volume: SYR

Response is of the following format:

space, space, f f f f f

IN d

Request to read the TTL logic level of the specified pin on the external 9 pin D-SUB connector. Valid pin numbers for input are:

6, 7, 8 and 9

If the pin specified is valid and if the pin level is high, response will be:

ON

If the pin level is low, response will be:

OFF

If the pin specified is invalid:

An 'Out Of Range' response will be given

OUT d = <ON|OFF>

Request to set the TTL logic level at the specified pin on the external 9 pin D-SUB connector.

Valid pin number for output is: 4

If the pin specified is invalid:

An 'Out Of Range' response will be given

Example:

Set pin 4 high:

OUT 4 = ON

Set pin 4 low:

OUT 4 = OFF

Pump Chain Commands

SEQ [<integer>] [<entry>]

Request to set or query programming sequences. <integer> is sequence number.

Default is Sequence 1. Command only applicable while pump is stopped.

Valid

sequence numbers, 'n', are 1 to 10. 'n' defaults to 1 wherever it is optional.

See program

examples.

Query entire program: SEQ

Example response:

```
SEQ 1: DISPENSE
75.000 ml/mn
43.155 ml
0:00:01 INTERVAL
  3 REPEAT
INFUSE
SEQ 2: PROFILE
100.00 ml/mn
150.00 ml
REFILL
SEQ 3: RESTART
```

Query program sequence [n]: SEQ n

Example response to the command 'SEQ 2' with the previous example's program:

```
SEQ 2: PROFILE
100.00 ml/mn
150.00 ml
REFILL
```

Query program sequence n's mode: SEQ [n] MOD

Response will be according to the following table:

<u>Response</u>	<u>Description</u>	<u>Response</u>	<u>Description</u>
STP	stop	RST	restart
PRO	profile	GOT	go to
INC	increment	EVN	event
DEC	decrement	PMP	pump
DIS	dispense	OUT	TTL out
PAS	pause		

Pump Chain Commands

Query data item of program sequence n:

<u>Command</u>	<u>Description</u>
SEQ [n] RAT	Query rate Response: f f f f f units Where units is one of the following: ml/mn ul/mn ml/hr ul/hr
SEQ [n] GOT	Query go to sequence number Response: <n>
SEQ [n] TGT	Query target volume Response: <float>
SEQ [n] INT	Query time interval Response: <time>
SEQ [n] RPT	Query repetition count Response: <float>
SEQ [n] OUT	Query output pin level setting Response: <ON/OFF>
SEQ [n] DIR	Query pumping direction Possible responses: INFUSE REFILL

Set mode of program sequence n:

SEQ [n] MOD mode
Where mode is as follows:

<u>Mode</u>	<u>Description</u>	<u>Mode</u>	<u>Description</u>
STP	stop	RST	restart
PRO	profile	EVN	event
INC	increment	GOT	go to
DEC	decrement	OUT	set output pin
DIS	dispense	PMP	pump
PAS	pause		

Set data item of program sequence n:

Set sequence's rate:
SEQ [<n>] RAT <float> [<units>]
Rate is of format: f f f f f

<u>Units are one of :</u>		<u>Description</u>
UM		µl/mn
UH		µl/hr
MM		ml/mn
MH		ml/hr

Pump Chain Commands

Set sequence's go to sequence number

SEC [<n>] GOT <n>

Set sequence's target volume:

SEQ [<n>] TGT <float>

Set sequence's time or target:

SEQ [<n>] INT <time>

Set sequence's repetition number:

SEQ [<n>] RPT <integer>

Set sequence's pumping direction:

SEQ [<n>] DIR <INF|REF>

Set sequence's output pin level

SEQ [<n>] OUT <ON|OFF>

VER Request for version of pumps embedded software.

Pump Chain Error Messages

Error messages are in the format:

<lf>, space, space, <message>, <cr>.

Where <message> is one of the following:

? Syntax error in a received command

NA Command not applicable at this time

OOR Control data is out of the operating range of the pump

Note: If assistance is needed in programming pump control via RS-232, call Harvard Apparatus Customer Service Support at 1-800-272-2775.

Appendix A: Syringe Diameters in mm

<i>Stainless Steel</i>		<i>Terumo</i>		<i>SGE Scientific Glass Engineering</i>	
<u>Size</u>	<u>Diameter</u>	<u>Size</u>	<u>Diameter</u>	<u>Size</u>	<u>Diameter</u>
8 cc	9.525 mm	3 cc	8.95 mm	25 µl	0.73 mm
20	19.130	5	13.00	50	1.03
50	28.600	10	15.80	100	1.46
100	34.900	20	20.15	250	2.30
200	44.755	30	23.10	500	3.26
<i>Becton Dickinson Plastic 'Plasti-pak'</i>		<i>Sherwood-Monoject Plastic</i>		1.0 ml 4.61 mm	
<u>Size</u>	<u>Diameter</u>	<u>Size</u>	<u>Diameter</u>	2.5 7.28	
1 cc	4.78 mm	1 cc	4.65 mm	5 10.30	
3	8.66	3	8.94	10 14.57	
5	12.06	6	12.70		
10	14.50	12	15.90		
20	19.13	20	20.40		
30	21.70	35	23.80		
50/60	26.70	60	26.60		
<i>Air-Tite 'All Plastic'</i>		<i>Popper & Sons, Inc. 'Perfektum' Glass</i>		<i>Hamilton-Microliter Series Gastight</i>	
<u>Size</u>	<u>Diameter</u>	<u>Size</u>	<u>Diameter</u>	<u>Size</u>	<u>Diameter</u>
2.5 cc	9.60 mm	0.25 cc	3.45 mm	0.5 µl	0.103 mm
5	12.45	0.5	3.45	1	0.1457
10	15.90	1	4.50	2	0.206
20	20.05	2	8.92	5	0.3257
30	22.50	3	8.99	10	0.460
50	29.00	5	11.70	25	0.729
<i>Unimetrics Series 4000 & 5000</i>		<i>Renfac</i>		50 1.031	
<u>Size</u>	<u>Diameter</u>	<u>Size</u>	<u>Diameter</u>	100 1.46	
10 µl	0.460 mm	2 cc	9.12 mm	250 2.3	
25	0.729	5	12.34	500 3.26	
50	1.031	10	14.55	1.0 ml 4.61 mm	
100	1.460	20	19.86	2.5 7.28	
250	2.300	30	23.20	5 10.3	
500	3.260	50	27.60	10 14.57	
1000	4.610			25 23.0	
				50 32.6	
				<i>Becton Dickinson Glass-All Types</i>	
				<u>Size</u>	<u>Diameter</u>
				0.5 µl	4.64 mm
				1	4.64
				2.5	8.66
				5	11.86
				10	14.34
				20	19.13
				30	22.70
				50	28.60
				100	34.90

Appendix B: Stainless Steel Syringes



Harvard High Pressure Stainless Steel Syringes

Harvard now offers a completely new line of Stainless Steel Syringes intended for high pressure applications with good resistance to most aggressive liquids. Wetted parts are #316 stainless steel or Viton. Syringes are available in 20, 50, 100 and 200 ml sizes with removable replaceable tips. Genuine SWAGelok® syringe to tube fittings are available in 1/16, 1/8 and 1/4 inch sizes. A luer lock end fitting is also available. All tips are interchangeable with all syringes (20 to 200 ml) in the series.

Both syringe barrel end plungers are #316 stainless steel. A Viton O-ring seal between top and end of the barrel insures against leakage. Syringes are guaranteed to be leak free for pressures up to 700 psi.

All syringes are supplied with inside diameter dimensions for use with Harvard microprocessor controlled pumps and rate charts for use with older 'classic' pumps. Replacement Viton O-Rings are available, as are the more chemically resistant Chemraz® O-Rings.

Harvard High Pressure 8 ml Stainless Steel Syringe

This syringe has been designed to utilize the high forces available in our syringe pumps to produce pressures up to 1500 psi. The syringe is constructed entirely of #316 stainless steel with two Chemraz® O-Ring seals, two Teflon O-Ring seals and 1.6 mm (1/16 in) SWAGelok® end. When used in the 70-2200 or 70-2201 Harvard PHD 4400, pressures in excess of 1500 psi are achieved for chromatography and process control applications.

Specifications

Volume	8 ml
Dimensions:	
Length of Barrel	16.5 cm (6-1/2 in)
Plunger Excursion	11.4 cm (4-1/2 in)
OD	1.6 cm (5/8 in)
ID	0.95 cm (3/8 in)
Max Test Pressure	4000 p.s.i
Working Pressure	1500 p.s.i

Cat. No. Product

55-3369 Harvard High Pressure 8 ml Stainless Steel Syringe

Replacement Parts

- 5013-089 Chemraz® O-Ring 20 ml
- 5013-090 Chemraz® O-Ring 50 ml
- 5013-091 Chemraz® O-Ring 100 ml
- 5013-092 Chemraz® O-Ring 200 ml
- 5013-087 Chemraz® Tip Seal O-Ring, all sizes
- 72-2472 Replacement Viton O-Ring 20 ml, 10
- 72-2473 Replacement Viton O-Ring 50 ml, 10
- 72-2474 Replacement Viton O-Ring 100 ml, 10
- 72-2475 Replacement Viton O-Ring 200 ml, 10
- 72-2476 Replacement Viton Tip Seal O-Ring, 10
- 72-2477 Stainless Steel Plunger Button to Adapt Syringe for Use with PHD 2000 HVP with Triple S Clamp

Harvard High Pressure Stainless Steel Syringes

Syringe Size	With SWAGelok® Diameter		With Luer Lock	
	1/16 inch	1/8 inch	1/4 inch	
20 ml	55-4199	55-4205	55-4209	55-4214
50 ml	55-4200	55-4206	55-4210	55-4215
100 ml	55-4201	55-4207	55-4211	55-4216
200 ml	55-4202	55-4208	55-4213	55-4217
Replacement Tips, Furnished with Sealing O-Ring				
	55-4220	55-4221	55-4222	55-4223

Appendix C: Min/Max Flow Rates

Minimum flow rates are taken from the smallest inside diameters and maximum flow rates are taken from the largest inside diameters of the syringes supplied by the eleven most widely used syringe manufacturers.

Nominal Minimum/Maximum Flow Rates for Various Syringes (Actual Limits will vary depending on manufacturer)								
Syringe Size	$\mu\text{l/hr}$		$\mu\text{l/min}$		ml/hr		ml/min	
	Min.	Max.	Min.	Max.	Min.	Max.	Min.	Max.
0.5 μl	0.0001	95.330						
1 μl	0.0002	190.70						
2 μl	0.0004	381.30						
5 μl	0.0010	953.17						
10 μl	0.0019					1.9013		
25 μl	0.0046					4.7752		
50 μl	0.0092					9.5511		
100 μl	0.0183					19.153		
250 μl	0.0454					47.532		
500 μl	0.0911					95.492		
1000 μl			0.0031			190.95		
1 ml			0.0033			205.30		
2 ml			0.0119			747.35		
2.5 ml			0.0076			476.21		
3 ml			0.010					11.231
5 ml			0.0208					21.781
10 ml			0.0301					31.486
20 ml			0.0523					54.804
30 ml			0.0673					70.518
50 ml			0.1019					106.76
100 ml			0.1740					182.40
140 ml			0.2106					220.82

Stainless Steel Syringes

Only Harvard's stainless steel syringes can be used with this pump. They can withstand high pressure applications and provide years of service.

Flow Rates for PHD 2000 HVP (High Volume and Pressure)

Rates listed are for single syringe

Syringe Size	Minimum	Maximum
20 ml	1.5 $\mu\text{l/hr}$	20 ml/min
50 ml	3.4 $\mu\text{l/hr}$	46 ml/min
100 ml	5.0 $\mu\text{l/hr}$	68 ml/min
200 ml	8.2 $\mu\text{l/hr}$	112 ml/min

Appendix D: PHD 2000 to PC Connection

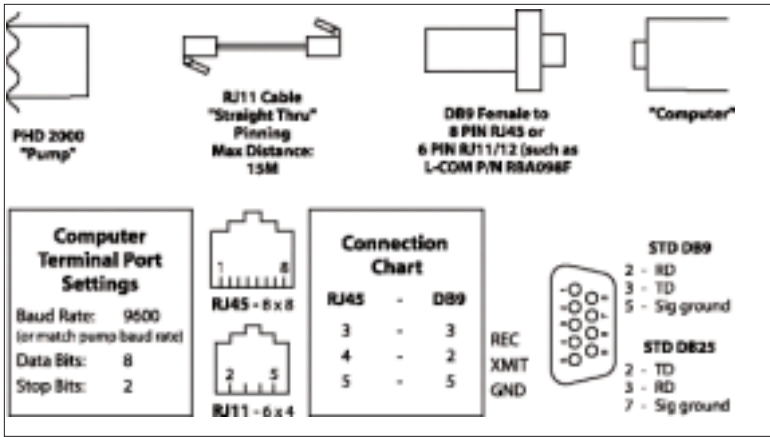


Figure 13. PHD 2000 to PC Connection

Appendix E: TTL Connector Specs

TTL Connections are via a 9 pin D-sub type connector. All inputs are internally pulled up to a logic high.

Pin Assignments

1. Vss (Ground Reference, logic low)
Falling edge starts/stops pump
2. Valve Control Output:
Low - Infusing
High - Refilling
3. Running Output Indicator
Low - Pumped stopped
High - Pump running
4. Programmable Output
5. Logic High Reference (5V, 400 μ A)
6. Foot Switch Input
7. Timer Input
Rising edge starts pump
Falling edge stops pump
8. Directional Control Input
Rising edge sets pump to infuse

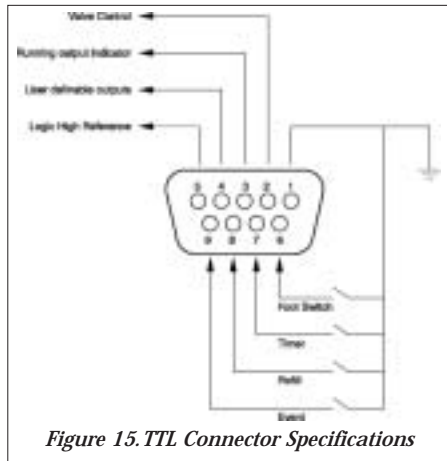


Figure 15. TTL Connector Specifications

- Falling edge sets pump to refill
9. Event Input
Falling edge triggers program event
- Logic Low:** 0V - 0.5V
Max. 2 mA current sink
- Logic High:** 2V - 5V
Max. 400 μ A current source

Appendix F: RS-232 Specifications

Pump Chain

- Pump Port 1 – Computer control side
- Pump Port 2 – Connection for remainder of pump chain
- Baud Rate – 1200, 2400, 9600 or 19,200
- Word Size – 8
- Parity – none
- Stop Bits – 2

Scale

Connect scale to Pump Port 1

Supported scales:

Mettler

- Baud Rate – 2400
- Word Size – 7
- Parity – even
- Stop Bits – 1
- Pause – 0

Sartorius

- Baud Rate – 2400
- Word Size – 7
- Parity – even
- Stop Bits – 1
- Code Settings:
 - 311 – Weigh only (models with PLUS performance package)
 - 211 – External print command without stability
 - 225 – 2400 baud
 - 234 – Even Parity
 - 432 – Beeper off

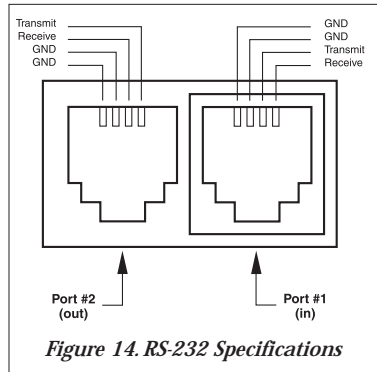


Figure 14. RS-232 Specifications

Ohaus

- Baud Rate – 2400
- Word Size – 8
- Parity – even
- Stop Bits – 1
- Auto print feature – off
- Stable data only – off
- Serial data frame selection – Fr.7

Printer

Serial printer with print buffer:

- Baud Rate – 2400
- Word Size – 8
- Parity – none
- Stop Bits – 2

Maintenance

PHD 2000 series pumps require no special maintenance other than keeping them clean by avoiding accidental spills of pumped material.

The two guide rods and the lead screw should be sparingly lubricated periodically. The guide rods and the lead screw should be lubricated with Magnalube-G R grease (Catalog #0606-086) provided with the pump.

Solvents of any type should never be used to clean the pump. A mild detergent solution may be used to clean the keypad.

Troubleshooting

Infusion Accuracy

To ensure infusion accuracy always use new syringe(s) and measure syringe bore diameter and enter actual dimensions in millimeters (mm). Additionally make sure that the guide rods and lead screw are properly lubricated.

RS-232 Difficulties

Verify that the baud rates and data framing parameters on all devices are the same. With the pump chain, a "Communication timeout" error on a computer is usually caused by the computer errantly handshaking on the RTS, CTS and DSR lines. Verify pins 4, 5 and 6, on the 25 pin connector, are jumped on the computer side of the cable.

PHD 2000 Display Messages

Pump Stalled

This indicates that the pusher block travel has been impeded. This may be caused by the syringe plunger hitting bottom, a kink in the tubing, syringe plunger binding or any situation requiring more force to the head of the syringe than the pump is capable of delivering. To restart the pump, remove the cause of stalling and press RUN. The pump will continue the interrupted procedure where it stopped. Also note, a large increase in the pumping rate could also stall the motor.

Out of Range

A value was entered or encountered in a pump program that was beyond the pump's limits. Also, verify that the Infuse Rate is not 0. Note that rates must be less than 42949. Express larger rates using different units.

EEPROM Data Lost

The values stored in the non-volatile memory have been lost. This should never occur under normal operation.

Replace NOVRAM

The non-volatile memory has failed. This should never occur under normal operation.

Appendix H: Accessories & Spare Parts

Auto Fill Accessory

To operate the Auto Fill feature, either the low pressure or high pressure valve assembly is required. Each assembly contains the solenoid valve, connector cable, power cable, and tubing and fittings to connect to the syringe. Operates at 115/230 VAC, 50/60 Hz via a selector switch.

Catalog No.	Description
Accessories	
55-7760	Daisy-Chain Cable
70-2022	Daisy-Chain Connector, 9 PIN
55-4145	Daisy-Chain Connector, 25 PIN
55-4144	Foot Switch
55-1146	'Auto Fill' Valve Box—Regular Pressure – to 25 psi
55-1145	'Auto Fill' Valve Box—High Pressure – 200 psi
59-7357	Manual Liquid Switch
59-7358	Automatic Liquid Switch—115 VAC, 60 Hz
59-7359	Automatic Liquid Switch—230 VAC, 50 Hz
59-7356	Replacement Tubing Connection Kit
59-7355	Replacement Reducing Adapters
61-0270	Continuous Flow Tube Segment**
72-0199	Remote Extension Cable, 1.5 m (5 ft)*
72-1405	Remote Ext. Cable, 9.1 m (30 ft)*
70-2022	PHD 2000 RS-232 Adapter, 9-pin to standard phone jack*
55-4202	SS Syringe, 200 ml, with 1/16 inch Swagelok®*
55-4208	SS Syringe, 200 ml, with 1/8 inch SWAGELOK®*
55-4213	SS Syringe, 200 ml, with 1/4 inch SWAGELOK®*
55-4217	SS Syringe, 200 ml, with Luer Lock*

* For PHD 2000 HVP Remote Syringe Pump

** For PHD 2000 Push/Pull Syringe Pump

Spare Parts

0606-086	Magnalube Lead Screw Lubricant
2400-243	Syringe Clamp
5012-005	3/32 Hex Key
5091-030	Syringe Retainer Knob
5100-055	T Knob - Withdraw and Anti-Siphon Bracket
5153-209	Power Cord
5400-001	PHD2000 Operators Manual

Appendix I: Kits and Upgrades

Catalog No.	Description
Kits	
70-2012	PHD 2000 6/10 Multitrack Upgrade Kit
70-2013	PHD 2000 4 x 140ml Multitrack Upgrade Kit
70-2014	PHD 2000 Microliter Rack Upgrade Kit
70-2015	PHD 2000 Anti-Siphon Kit (for Infusion Pump)
Upgrades	
70-2016	PHD 2000 Infusion to Infuse/Withdraw*
70-2017	PHD 2000 Infuse/Withdraw to Programmable*
70-2018	PHD 2000 Infusion to Programmable*

**Factory Installed Only*

Appendix J: Custom Applications

The Harvard PHD 2000 Syringe Pumps lend themselves to a multitude of OEM industrial applications. For all types of custom pumping or pilot plant applications, the pump's programming capabilities can be enhanced. Please contact the Harvard Development Group if we can be of help.

Appendix K: Symphony

Symphony, the Windows[®] Pump Manager software, allows you to control your PHD 2000 pump from your remote computer. With Symphony, you can orchestrate all pump functions from a virtual control panel via the serial communications link built into every PHD 2000 pump.

Catalog Number	Product
-----------------------	----------------

70-3000	Symphony — The Windows™ Pump Manager
---------	--------------------------------------

Harvard Apparatus pumps are distributed
on a non-exclusive basis by:

INSTECH

Instech Laboratories, Inc.

Headquarters:
5209 Militia Hill Road
Plymouth Meeting, PA 19462
TL (800) 443-4227
TL (610) 941-0132
FX (610) 941-0134

European sales office:
P.O. Box 2
9950 AA Winsum
The Netherlands
TL 0031-(0)594-696123

www.instechlabs.com/Pumps/syringe/