



## Circulator Set-up



Insulated tubing connections to circulator.



Align thermal block for connection.




Insert fluid Quick Disconnect to thermal block.



Set desired circulator temperature setting.

## Insert and Thermal Block Set-up

The following insert and thermal block photo sequence details how the plate fits the insert and then placement of both onto thermal block.

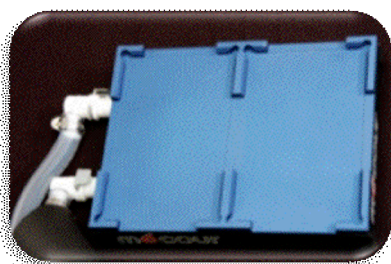

Two-plate "Open Well" thermalblock.



PCR Plate & thermal insert.



***Insert and Thermal Block Set-up continued***



Locate plate on PCR thermal insert.



Load plate/insert on thermal block.




Easy loading of reagent reservoir to Thermal Insert




Flexible changing of consumables



Easy access and loading

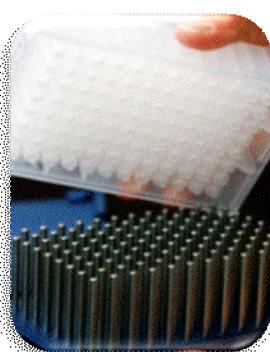

Deep well plate

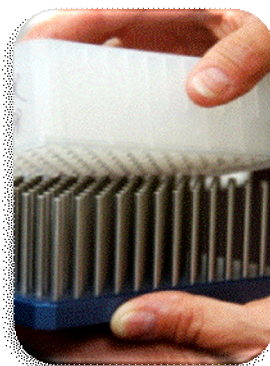

Loading DWP to thermal insert



Ready for Thermal Block placement