



DISCOVERY

L I F E S C I E N C E S

Science at your Service.™

Lung Dissociated Tumor Cell Thawing and Culturing Procedure

Safety:

- Primary tumor cells should be contained in a Class II biological safety cabinet and handled using Biosafety Level 2 (BSL-2) work practices and facilities. (2)

Cell Culture Media and Antibiotics:

- **Recommended cell culture media (3):**
 - DMEM-F12 medium with L-glutamine
 - 5 mM HEPES
 - 0.1% sodium bicarbonate
 - 0.4% BSA or 10% FBS
- **Antibiotics:**
 - Normocin (Invivogen; 100 µg/mL)
- **Supplements:**
 - 20 ng/ml EGF
 - 10 ng/ml bFGF
 - Supplements recommended for unthawing samples
 - DNase I (1 mg/mL), resuspended in PBS without calcium or magnesium and filter-sterilized

Unthawing Procedure:

1. On the day of the thaw, pre-warm 19 mLs of recommended cell culture media in a 37°C incubator.
2. Divide the media into two 15-mL conical tubes (1 tube with 9 mLs of media; 1 tube with 10 mLs of media).
3. Quickly thaw the DTC sample in a 37-40°C water bath until 2mm crystals remain. Move sample through the water to speed thawing.
4. Slowly add the 1-mL DTC sample to 9 mLs of pre-warmed media. Gently mix by inversion (do not vortex).

Lung Dissociated Tumor Cell Thawing and Culturing Procedure - Continued

5. Centrifuge the conical tube at 300 x g for 5-10 minutes (no brake). Carefully remove the supernatant.
6. After the initial wash to remove DMSO from the sample, gently resuspend the pellet in the remaining 10 mLs of pre-warmed cell culture media.
7. If the cells clump together, perform the following:
 - a. Add DNase I (1 mg/mL) to the cell culture media.
 - b. Incubate at 37°C for 5-10 minutes to facilitate digestion of released DNA.
 - c. Centrifuge the conical tube at 300 x g for 5 - 10 minutes (no brake).
 - d. Carefully remove the supernatant containing DNase I.
 - e. Gently suspend the pellet in 10 mLs of pre-warmed cell culture media. Go to the next step.
8. Allow cells to rest for at least 1 hour in the media at 37°C before checking the viability and plating.
9. Check the viability and cell number of the sample.
10. Plate the cells according to assay requirements.

Culturing Primary Cells:

1. For sphere formation assays, using untreated tissue culture flasks is recommended to reduce cell adherence (4).
2. The media recipe listed above is a base media for thawing and acclimating lung tumor cells, in preparation for cell culture.
3. Additional supplementation will be required to expand lung tumor primary cells in vitro (3), depending on the end-goal of the assay. Please see the literature for supplementation recommendations, depending on the specific needs of the cell types targeted for expansion.

References:

1. Chan, L. L.; D. J. Lavery; T. Smith; P. Nejad; H. Hei; R. Gandhi; D. Kuksin; J. Qiu. 2013. Accurate measurement of peripheral blood mononuclear cell concentration using image cytometry to eliminate RBC-induced counting error. *J Immunol Methods*. 388:25-32.
2. Centers for Disease Control and Prevention. 2009. Biosafety in Microbiological and Biomedical Laboratories (BMBL). U.S. Dept. of Health and Human Services. 5th ed.
3. Eramo, A.; F. Lotti; G. Sette; E. Pilozzi; M. Biffoni; A. Di Virgilio; C. Conticello; L. Ruco; C. Peschle; R. De Maria. 2008. Identification and expansion of the tumorigenic lung cancer stem cell population. *Cell Death and Differentiation*. 15:504- 514.
4. Pan, J.; Q. Zhang; Y. Wang; M. You. 2010. 26S proteasome activity is down-regulated in lung cancer stem-like cells propagated in vitro. *PLoS One*. 5:e13298