

SCIENCE. LOGISTICS. CERTAINTY

High Volume Shipment: Loading

Sample Provider

How to Load Non– Infectious Samples Page 1





SCIENCE. LOGISTICS. CERTAINTY.

High Volume Shipment: Loading

Sample Provider

How to Load Non– Infectious Samples Page 2

Section D	F	Packaging the Shipment
Step D1 Replace the shippe	r plug. Remove the Leg 2 ship- ping pouch from docu- ment protec- tor attached to the dewar handle in- side of the container.	Step D3 Step D3 Remove any classification labels from the shipping pouch and place on the metal diamond. Image: Step D3 Image: Step D3
		 Step D5 Once all documents are loaded: 1. Close the shipping pouch and remove sticker backing. 2. Place shipping pouch on metal plate. 3. Close lid.
Step D6 Image: Constraint of the state of th	Step D7Always secure Latches with Zip Ties prior to ship- mentSee these instructions for Zip Tie and Tamper Proof Tag procedures9210 CXHVO Latch Zip Tie Procedure9211 CXHVO Tamper Proof Tag Procedure	For those shipments containing regulated dangerous goods/hazardous materials, the shipper is responsible for correctly preparing the shipment according to the current International Air Transport Association (IATA) and International Civil Aviation Organization (ICAO) dangerous goods regulations. This includes correct identification, classification, packaging, dangerous goods markings and labeling as well as completion of all pertinent and required documentation. The shipper is the customer who is preparing the shipment at each leg in the transportation of the dry dewar packaging.

24-HOUR HELP LINE (949) 470-2305

CryoPort, Inc. • 17305 Daimler St. • Irvine, CA 92614 Phone: 1-949-470-2300 Fax: 1-949-470-2306 www.cryoport.com • E-mail Contact: cs@cryoport.com