

Sample Provider

How to Load Infectious
Samples
Page 1

High Volume Shipment: Loading the SafePakXL™

Section A

Step A1



Once the dewar has arrived, remove old shipping pouch.

Receiving Shipper

Step A2



Unlock latch by turning the key to the left.

The latch will "unhook," from the top of the dewar.






Open the lid to expose the dewar.

Section B

Preparing the SafepakXL™

(Please refer to Cryostrap[™] Instructions)

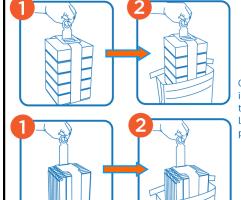

Open Bag and insert one rack into top opening.
Lower rack completely into bag.



X Detail A

Flatten top section of pouch. Remove cover from adhesive strip, fold along crease (Detail A) and seal bag.



Fold tabs around edges and seal.

Section C

Form: 9189 v2 Effective: 03/6/17

Loading Samples for Shipment

IMPORTANT NOTE: Commodity and accessories must be pre-conditioned prior to loading into the sample chamber. A 20% reduction of holding time by can occur if the materials are not preconditioned prior to loading into the shipper.

Load the SafepakXLTM into the Dewar



Allow it to seat in the bottom of the Dewar



Install the Vapor Plug and allow it to seat down completely



24-HOUR HELP LINE (949) 470-2305

Instructions Continued on Back of Sheet.



SCIENCE. LOGISTICS. CERTAINTY.

Sample Provider

How to Load Infectious
Samples
Page 2

High Volume Shipment: Loading

Section D

Packaging the Shipment

Step D1

Replace the shipper plug.



Step D2

Remove documents from the shipping pouch # 1 from the



Step D3

Remove any classification labels from the shipping pouch and place on the metal diamond.

Step D4

International Shipments ONLY

Remove the Commercial Invoices located in the shipping pouch.

- 1. Sign and Date.
- Return all documents being shipped behind the Air Waybill.

 Including any Parmits Forms
 - <u>Including any:</u> Permits, Forms, Licenses, etc.



Step D5

Once all documents are loaded:

- Close the shipping pouch and remove sticker backing.
- 2. Place shipping pouch on metal plate.
- Close lid.







Step D6



Lock latch by turning the key to the right.

Step D7



For those shipments containing regulated dangerous goods/hazardous materials, the shipper is responsible for correctly preparing the shipment according to the current International Air Transport Association (IATA) and International Civil Aviation Organization (ICAO) dangerous goods regulations. This includes correct identification, classification, packaging, dangerous goods markings and labeling as well as completion of all pertinent and required documentation. The shipper is the customer who is preparing the shipment at each leg in the transportation of the dry dewar packaging.

24-HOUR HELP LINE (949) 470-2305 CryoPort, Inc. • 17305 Daimler St. • Irvine CA 92614-5510 Phone: 1-949-470-2300 Option 1 Fax: 1-949-470-2306 www.cryoport.com • E-mail: cs@cryoport.com

Form: 9189 v2 Effective: 03/6/17