




#### Step 1

Open the SmartpakII™ Housing located on the Vapor Plug of the CXHVO shipper by unlatching the black latch and pivoting the lid upwards

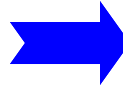

#### Step 2

Partially lift the SmartpakII™ from the foam cradle.  
Disconnect the thermocouple from the mini-USB port.

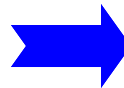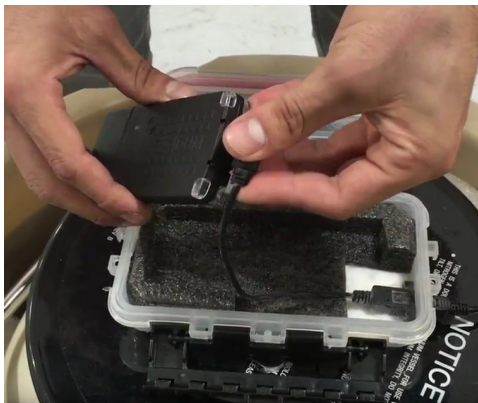

#### Step 3

Lift the SmartpakII™ up out of the container.




To Reassemble Unit follow this Procedure in Reverse Order.




**24-HOUR HELP LINE  
(949) 470-2305**

**Instructions Continued  
on Back of Sheet.**


#### Reset

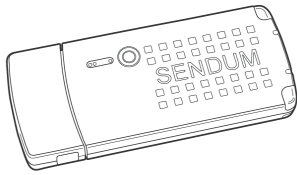


SmartpakII™ removed from packaging

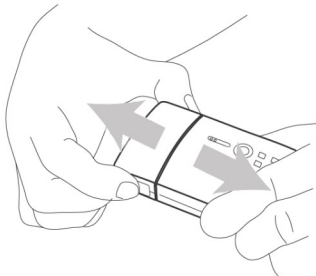

Push Buttons on both sides of the unit to unlock

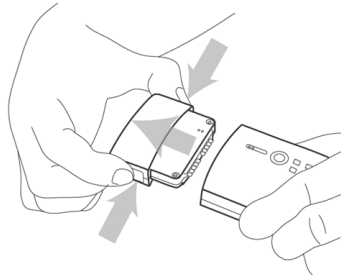


Pull sections apart and hold for a count of approx. 20 seconds

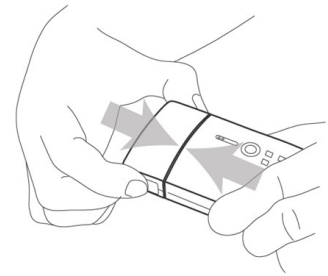


Push sections together until you feel and hear the sections click together


#### Replace

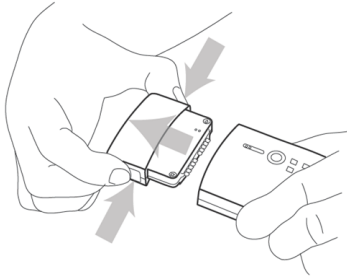

Detach Control Module and reattach it to a Fresh Battery Pack to the new battery pack

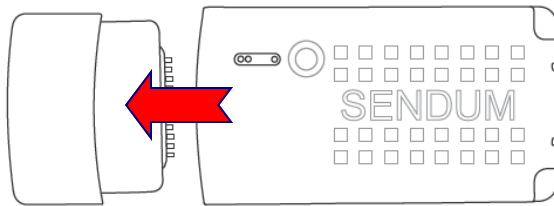


Depleted Battery


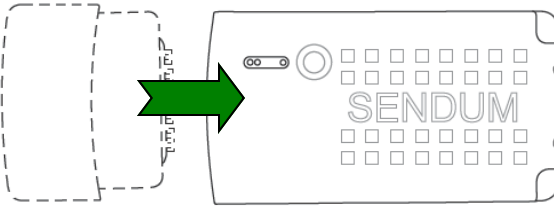


Fresh Battery

#### Recharge

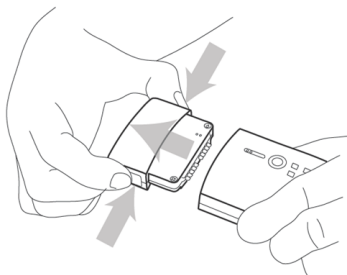

Detach Control Module and attach it to a Mini USB Charger for 6 hours.

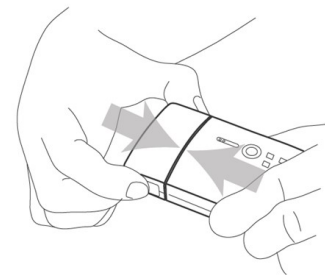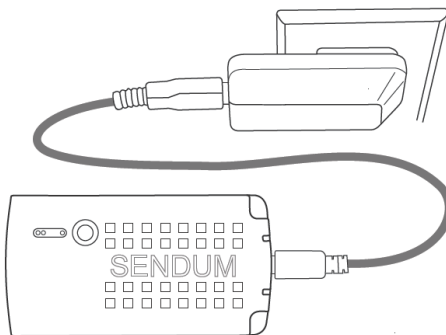

Reattach Control Module to Charged Battery Pack

**24-HOUR HELP LINE**  
**(949) 470-2305**

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614

Phone: 1-949-470-2300 Fax: 1-949-470-2306

www.cryoport.com • E-mail Contact: cs@cryoport.com