

PERSONAL PROTECTIVE EQUIPMENT:


- Cryogenic Gloves
- Safety Goggles

Step 1



Unlock latch by turning the key to the left.

The latch will release from the top of the dewar.

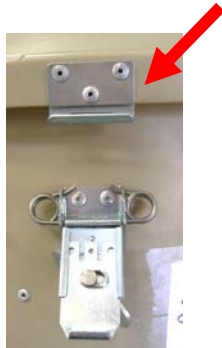

Open the lid to expose the dewar.

Step 2

Remove the vapor plug and rest it sideways, behind the dewar and against the brown top.




Step 3

1. Slowly pour Liquid Nitrogen into opening. (Cryoport Client Care will provide information on the amount of LN2 to use for your specific use case.)



Step 4

Replace the vapor plug and let the Dry Shipper stabilize, undisturbed, for at least 30 minutes, Before handing it over to carrier

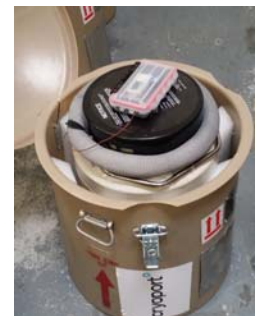

Step 5

Complete the In Field Recharging form and send it to cs@cryoport.com

cryoport In-Field Recharging Verification Form Rev. 4/2017

Re-Charging Date: _____
 Re-Charging Time: _____
 Location: _____
 Name of Person Recharging Dewar: _____
 Name of Company: _____
 Location of Dewar: _____

Pre-Charging Information
 Please indicate if the dewar was previously used for LN2 recharging:
 Dewar contains commodity Dewar does NOT contain commodity
 Dewar is COLD Dewar is WARM

Weight of the Shipper Prior to Recharging: _____

Recharging Information
 Amount of LN2 added to Dewar: _____

Post-Charging Information
 Please indicate if the dewar was previously used for LN2 recharging:
 Dewar contains commodity Dewar does NOT contain commodity
 Weight of the Shipper After Recharging: _____

Signature of Person Recharging Dewar: _____

**24-HOUR HELP LINE
(949) 470-2305**

CryoPort, Inc. • 17305 Daimler St. • Irvine, CA 92614

Phone: 1-949-470-2300 Fax: 1-949-470-2306

www.cryoport.com • E-mail Contact: cs@cryoport.com