

Standard Shipment: CXHVO Data Logger Access

Sample Receiver


How to Access LCD Data Logger—CXHVO

Step 1



Open Shipper.

Step 2a

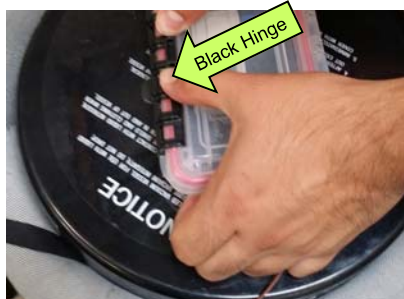

Unlatch Data Logger box using thumbs to press against the black hinge.

Step 2b



Properly unlatched Data Logger box.

Step 4



Open the Data Logger Box. If you are not downloading data, Go to Step 5a. If your Data Station Software is installed on a nearby computer, Go to Step 5a. Otherwise Go to Step 5b.

Step 5a




Remove the data logger from the box. You can now follow the Data Station instruction to download temperature data, if needed

Step 5b



If your Data Station Software is not installed on a nearby computer, disconnect Thermocouple from Data Logger. You can now follow the Data Station instruction to download temperature data.

Step 7

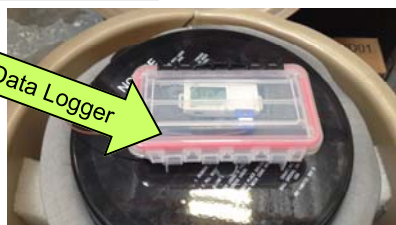

After downloading the data, return the data logger to the Data Logger Box. Re-attach the Thermocouple if required.

Step 8

Follow the Unloading Instructions to complete the receiving process and return the shipper to Cryoport.



**24-HOUR HELP LINE
(949) 470-2305**

CryoPort, Inc. • 17305 Daimler St. • Irvine, CA 92614
Phone: 1-949-470-2300 Fax: 1-949-470-2306
www.cryoport.com • E-mail Contact: cs@cryoport.com