


As Easy as 1 2 3

1 Receive

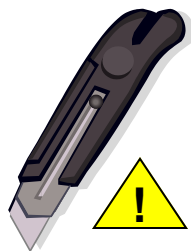

Use caution when using blade to open brown carton. Do not cut through box.

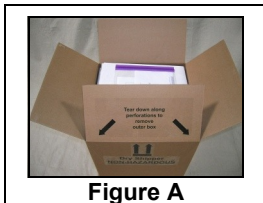


Figure A



Figure B



Figure C

Remove and discard outer BROWN shipping carton by grasping the top brown carton flap marked with arrows (fig. A) and pulling down along perforations on side of brown carton (fig. B).

Brown carton should be discarded (fig. C). Do not use for any further shipping!

2 Prepare



UNLOAD & RE-LOAD SAMPLES


See "Unload Samples" instruction sheet for full details on the back of this sheet.

(Loading Instructions are located on Flap "A" of white carton)



Flap "A" has paperwork already in place to ship the samples.

3 Ship

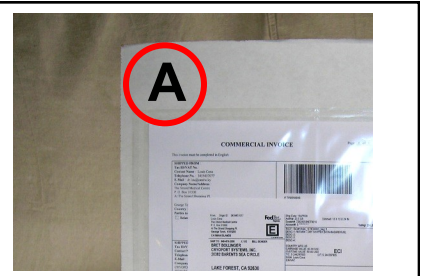

When preparing to ship, make sure flap "A" is on **TOP** of white carton! Be sure you followed "How to Load Samples" instruction sheet.

(Loading instructions are located on the back side of flap "A")



Seal top of white carton with packaging tape.

Carton is ready for pick-up.

24-HOUR HELP LINE
(949) 470-2305

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614
Phone: 1-949-470-2300 Fax: 1-949-470-2306
www.cryoport.com • E-mail: contact: cs@cryoport.com

Step 1




Wear Safety Gloves.
Hold shipper plug in one hand.
Hold blue strap with other hand.
Pull shipper plug upwards and out.

Step 2




Put shipper plug to one side.
Pull upwards on blue straps.
Remove sample bag completely
And place on a flat surface.

Step 3

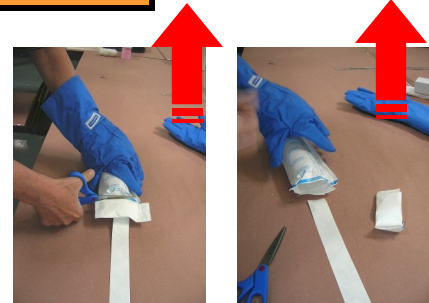


Hold bag with one gloved hand. Use
scissors to cut off end of Tyvek bag.
Remove end of bag completely to re-
veal sample container(s).

Step 4

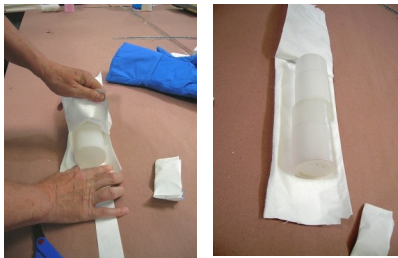

Tear backwards on Tyvek bag.
Remove bag completely.

Step 5

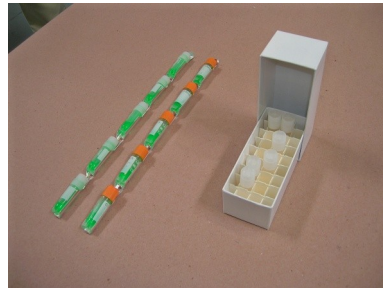

Retrieve sample container.

Step 6



Replace shipper plug.
Push plug fully into place.

***(Loading instructions are located on the back side of
flap "A")***

**24-HOUR HELP LINE
(949) 470-2305**

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614
Phone: 1-949-470-2300 Fax: 1-949-470-2306
www.cryoport.com • E-mail: contact: cs@cryoport.com