


STANDARD SHIPMENT: CXHVO LATCH TAMPER PROOF TAG PROCEDURE FOR THE DEWAR BODY


PRIMARY METHOD

STEP 1

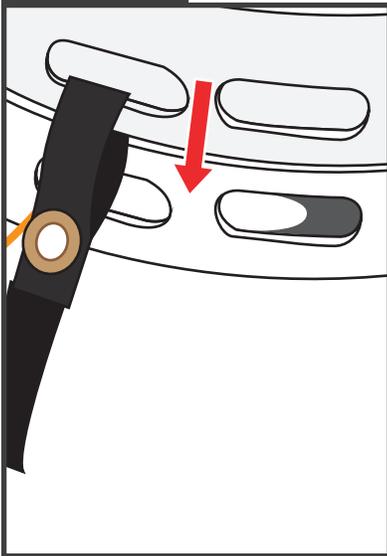


Align the slots in the vapor plug lid and the dewar. Close the lid.

STEP 2

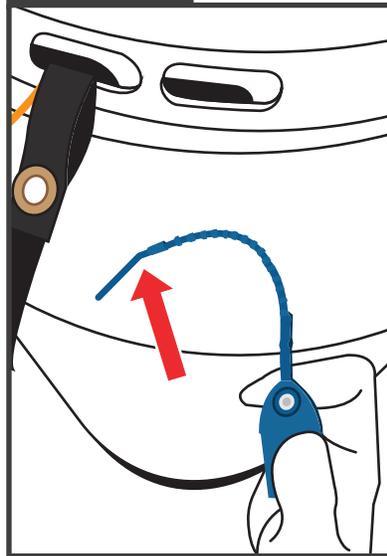


Take Tamper Proof Tag and slightly bend near the tip to aid in inserting the tag through the slots.

STEP 3

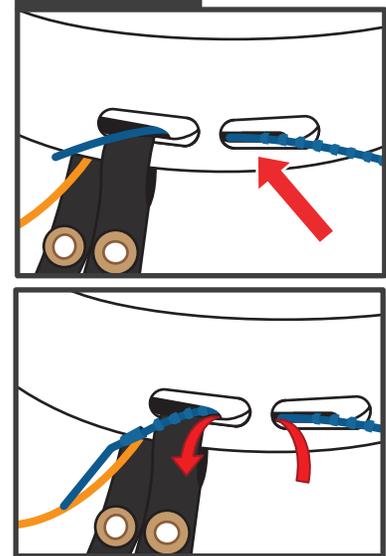

Insert Proof Tamper Tag through Vapor Plug Lid & Dewar slots.

STEP 4

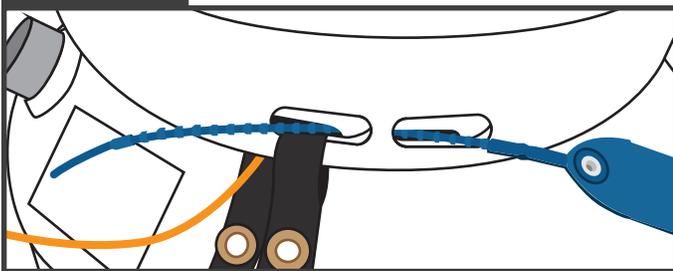


Pull the Tamper Proof Tag through slots until it is centered in the slots.

STEP 6

Install Zip Ties to secure enclosure latches per Instructions 9210.

STEP 5

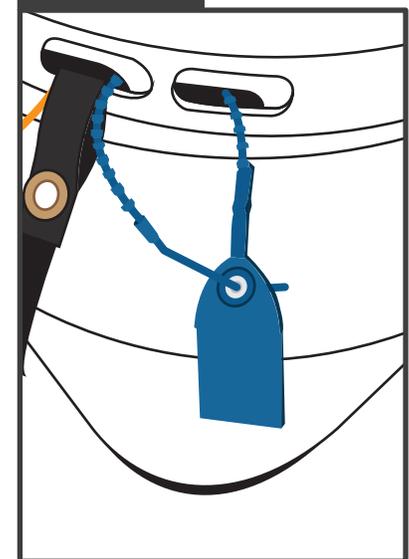


Insert Tamper Proof Tag tail through the Tamper Proof Tag's flag retention hole & pull tight.

IF THE DRY SHIPPER DOES NOT EMIT VAPOR WHEN PLUG IS REMOVED
PLEASE CONTACT CUSTOMER SERVICE IMMEDIATELY, BY CALLING (949) 470-2305 OR EMAIL CS@CRYOPORT.COM


STANDARD SHIPMENT: CXHVO LATCH TAMPER PROOF TAG PROCEDURE FOR THE SHIPPER ENCLOSURE

ALTERNATE METHOD

STEP 1

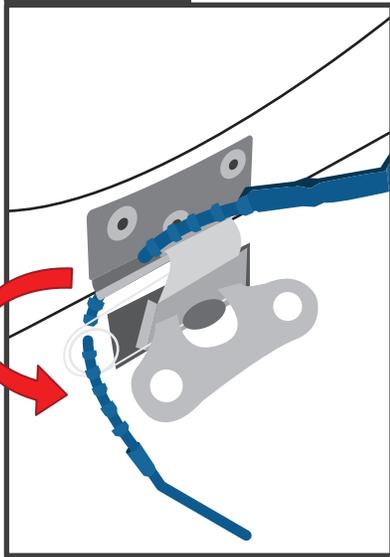


Insert Tamper Proof Tag down through Lid Latch Catch.

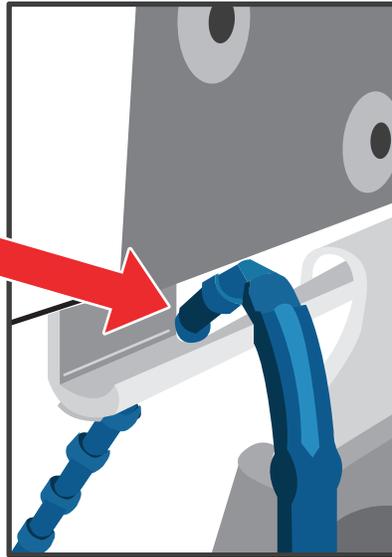


Assure that it is through the Latch Catch Slot.

STEP 2

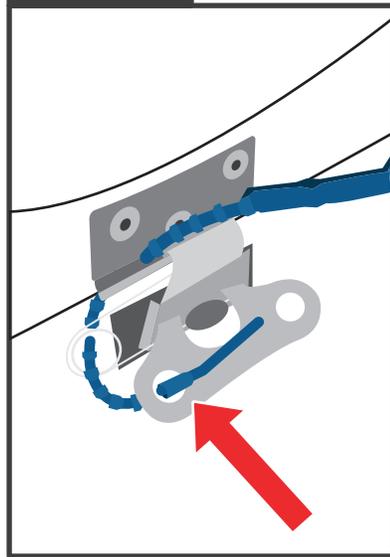

Thread Tamper Proof Tag Up through Body Latch Turnbuckle hole.

STEP 3

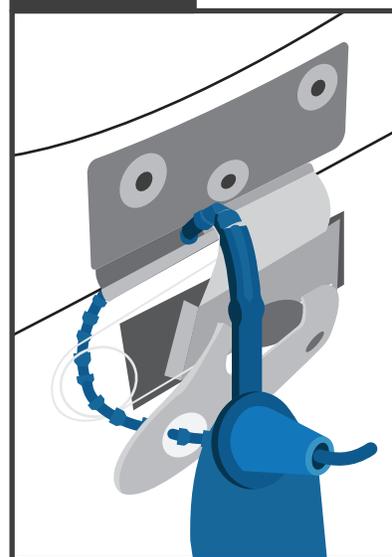

Insert Tamper Tag through Tamper Tag Flag retention hole & Pull Tight.

IF THE DRY SHIPPER DOES NOT EMIT VAPOR WHEN PLUG IS REMOVED
PLEASE CONTACT CUSTOMER SERVICE IMMEDIATELY, BY CALLING (949) 470-2305 OR EMAIL CS@CRYOPORT.COM

24-hour Help Line
(949) 470-2305

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614
Phone: 1-949-470-2300 | Fax: 1-949-470-2306
www.cryoport.com • e-mail contact: cs@cryoport.com