


How many specimens can TruFade accommodate?

The carousel inside TruFade accommodates 9 tri-sided specimen holders.

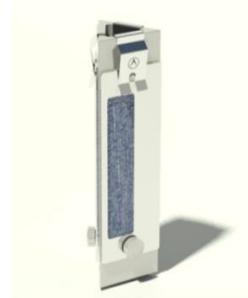

The holders can be fitted with an assortment of masks which allow the user to make comparisons between exposed and unexposed areas of the specimen and any control materials used:

Masks for ISO 105-B02: Method 1 are supplied with the [TruFade](#)



Single slot mask giving 2/3 cover



Double slot masks giving 1/3 cover

Masks for ISO 105-B02 : Method 2 are available as an optional accessory



Single slot mask giving quarter ($\frac{1}{4}$) cover



Single slot mask giving half ($\frac{1}{2}$) cover



Single slot mask giving three-quarters ($\frac{3}{4}$) cover

Each of the 9 holders have 3 sides that can accommodate a maximum of 10 specimens of ~1cm per side (30 specimens per holder) using the masks above.

TruFade can hold 270 specimens before blue wools or SolarSens™ are added.

Normally, SolarSens™ will occupy one of the spaces.

James Heal can also supply masks for compliance with AATCC methods.

These require larger specimens and can be fitted on any of the 27 sides, some of which may be replaced by blue wools or SolarSens™.

Special AATCC masks are available to ensure that the exposed areas conform to AATCC specifications.



Double slot mask with exposed areas
30x30mm



Double slot mask with exposed areas
30x15mm