

Thermostat Check – AccuDry2 417

WARNING: Any maintenance work undertaken on a James Heal instrument must be done so by a trained and competent person. Where electrical work is detailed, this should only be completed by a qualified engineer. James Heal is not responsible for damage or injury arising from work carried out by any person not qualified to do so. If in doubt, please contact James Heal or your local James Heal Agent.

Disconnect the Accudry2 from the electrical supply.

Using a 7mm nut-driver or socket remove the 3 screws in the centre of plastic cover.

Lift off the plastic cover.



Remove the 7 screws holding the metal duct at the rear of the dryer.



Remove the 1 screw holding the heater assembly to the dryer.



You will see 2 thermostats at the back of heaters.

Check the electrical resistance of these thermostats.
Both should be zero Ohms.

If any measure open circuit then they have failed.

