## Hydrosense Multi Legionella Water Test

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
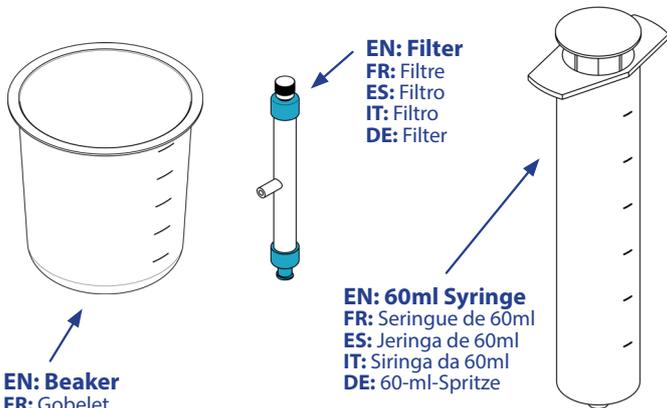
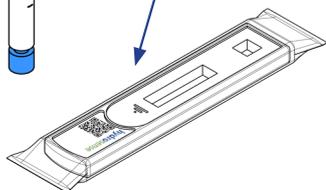
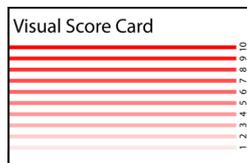
# 1

**EN: Collect 100ml of water from any water source using the beaker.**

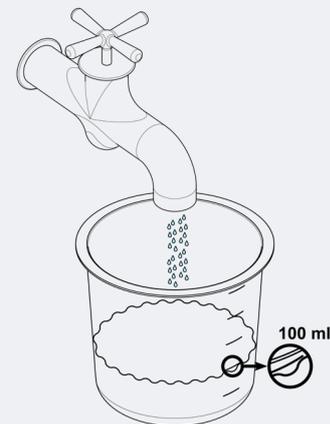
**FR: Prélevez 100ml d'eau de n'importe quel point d'eau à l'aide du gobelet.**  
**ES: Recoger 100ml de agua de cualquier ubicación en el vaso de precipitados.**  
**IT: Prendere 100ml di acqua da qualsiasi erogatore idrico utilizzando il becher.**  
**DE: Entnehmen Sie aus jeder Wasserquelle 100ml Wasser mit dem Becherglas.**

# 2

**EN: First, take the filter from its packaging and tighten the end cap.**

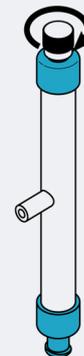**FR: Tout d'abord, sortez le filtre de son emballage et resserrez le capuchon.**  
**ES: En primer lugar, extraer el filtro de su embalaje y apretar el tapón.**  
**IT: Per prima cosa rimuovere il filtro dalla confezione e stringere bene il cappuccio all'estremità.**  
**DE: Nehmen Sie zuerst den Filter aus der Verpackung und ziehen Sie die Endkappe fest.**

# 4

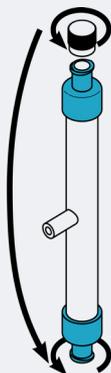
**EN:** Remove the end cap from the filter and place it on to the opposite end of the filter.

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# 5

**EN:** Take the 1ml recovery buffer syringe from the kit. Remove the blue cap.

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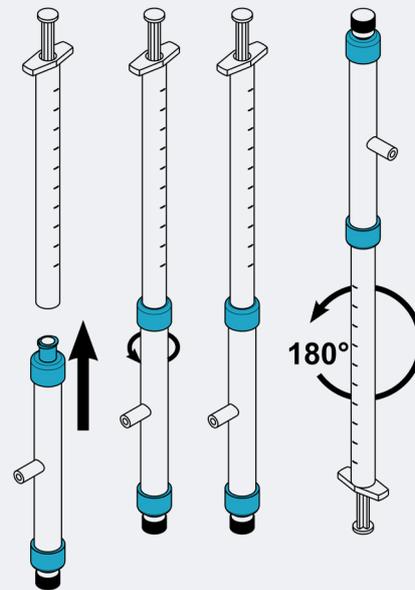
# 6

**EN:** Keeping the filter upright, attach the 1ml syringe to the open end of the filter. Flip the filter and 1ml syringe 180°.

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
# 7

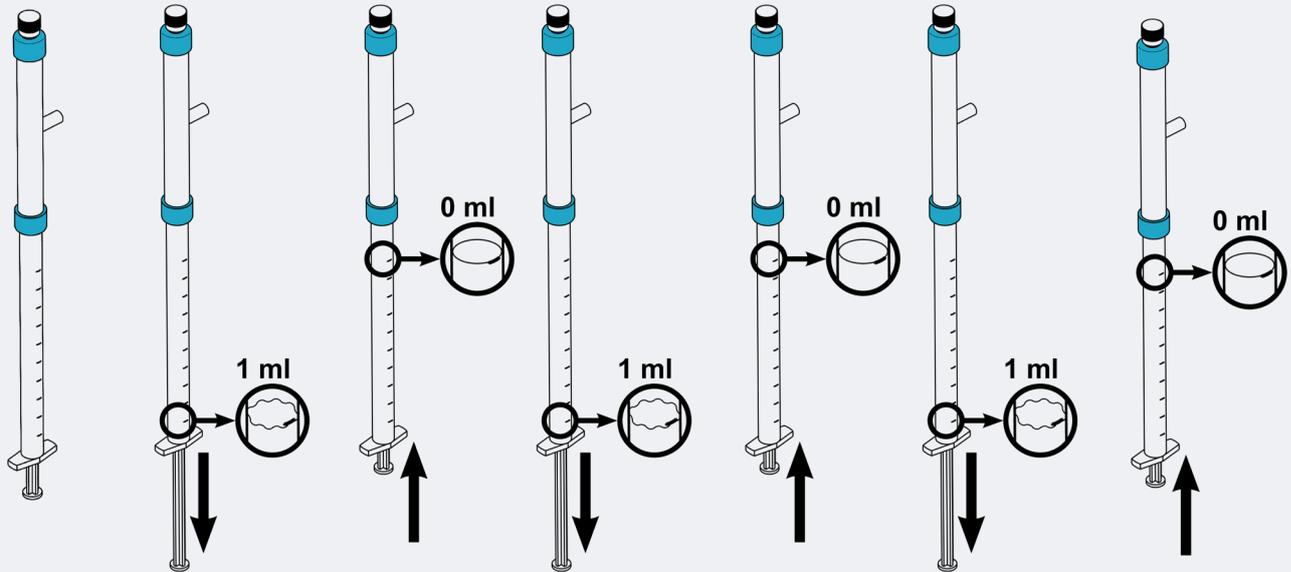**EN: Pull the syringe back to the 1ml mark. Push the liquid all the way back into the filter. Repeat this step 2 more times.**

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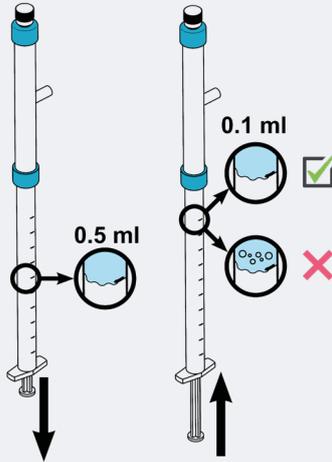
# 8

**EN: Pull the syringe back to the 0.5ml mark. Push the plunger in slowly to the 0.1ml mark to collect 0.1ml of sample exactly.**

*If there are any air bubbles in your sample, repeat the last step until there are no air bubbles in your 0.1ml sample.*

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# 9

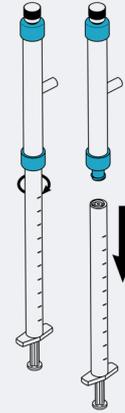
**EN: Detach the syringe from the filter, place the filter to one side.**

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# 10

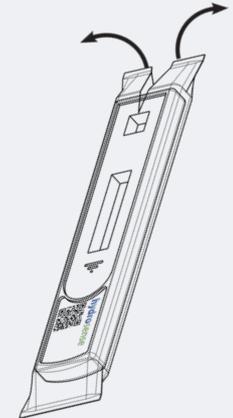
**EN: Take the test strip out of its foil packaging. Place the test on a flat surface.**

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# 11

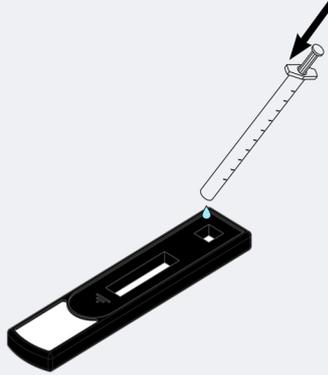
**EN: Gently add your 0.1ml of sample to the sample port on the test strip.**

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# 13

**EN: After 25 minutes (and before 30 minutes) read your Hydrosense Legionella test.**

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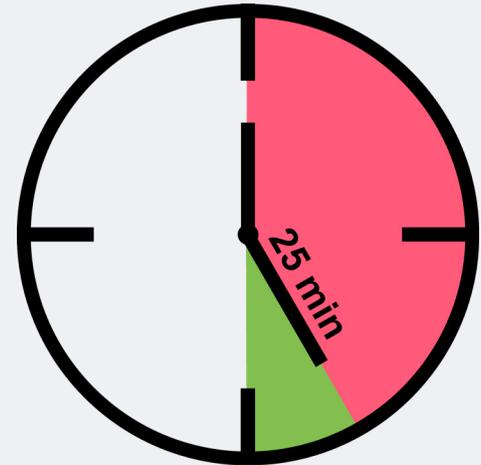
**EN: The test is composed of two lines; a test line (TL) and a control line (CL). If both lines are present, even if the test line is very faint, then *Legionella pneumophila* serogroup 1 has been detected above the cut-off of the test kit (>1,000 CFU/L).**

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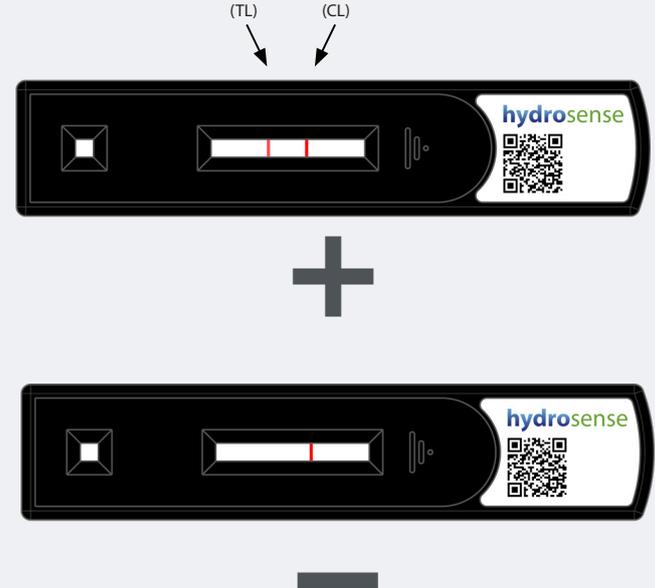
**EN: If no test line is present at the test line position but there is a control line, then the test is negative and *Legionella pneumophila* serogroup 1 has not been detected.**

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
**EN: Tests without a control line, or that look like any of these tests, are invalid.**

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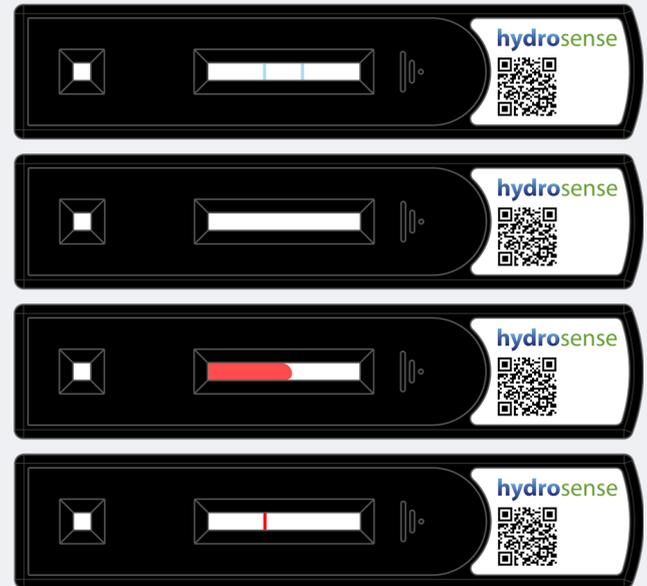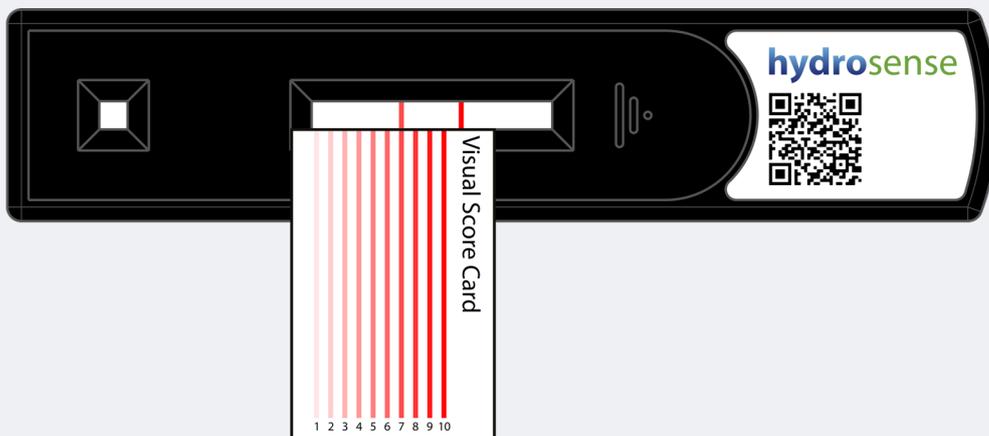
**EN: Read the Hydrosense test using the Hazard index card provided.**

**Use the visual score card to read the Hazard Index by matching the intensity of the test line on the test strip with the intensity of the lines on the visual score card.**

**A hazard index score of 1 or greater is equal to or greater than 1,000 CFU per litre and a hazard index score of 6 or greater is equal to or greater than 10,000 CFU per litre.**

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## THE SMALL PRINT

### EN: Limit of Detection:

1,000cfu/L *L. pneumophila* sg1 (when 100ml sample filtered). Filtration and test performance may be affected by suspended solids. Bacteria recovery post filtration can vary (<10-100%).

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### EN: Test Operating Limits:

- Use on samples between 10-45°C (50-113°F).
- Evaluated on samples that filter in <5 minutes.
- Do not use on systems treated with biguanide based or THPS biocides.

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### EN: Specificity:

- Non-reactive with wide range of environmental bacteria
- Weak cross reactions with other *L. pneumophila* sg's and *S. aureus* at high concentrations.

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### EN: Storage:

- Store at room temperature (18-22°C).
- Do not freeze.
- Do not use after expiry date.
- Do not use if test wrapper perforated.

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**EN: Disclaimer:**

Hydrosense Ltd makes no warranties or representations regarding performance of the products, or that the products are merchantable or fit for a particular purpose. Hydrosense expressly disclaims all other warranties and representations, express or implied, or which arise by operation of law or otherwise.

HYDROSENSE is a trademark of Hydrosense limited, used worldwide and registered in various territories.

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[©Hydrosense Ltd](http://Hydrosense Ltd)